



KERN & Sohn GmbH

Ziegelei 1
72336 Balingen-Frommern
Germany

+0049-[0]7433-9933-0
+0049-[0]7433-9933-149
info@kern-sohn.com

Installation instructions Tripods

KERN DS, CDS

Type
Version 1.0
2025-03
GB

TDS-A04-B
TDS-A05-B



TDS-A04-B_TDS-A05-B-IA-e-2510



Tripods
Version 1.0 2025-03
Installation instructions

Table of contents


1	Technical data.....	3
1.1	Dimensions.....	3
2	Scope of delivery	6
3	Installation	7
4	Appendix.....	10

1 Technical data




KERN	TDS-A04-B	TDS-A05-B
Tripod height (mm)	330	600

1.1 Dimensions




- Dimensions in mm
- Drawings on the next page
- Drawing of the adapter plate at the end of the document

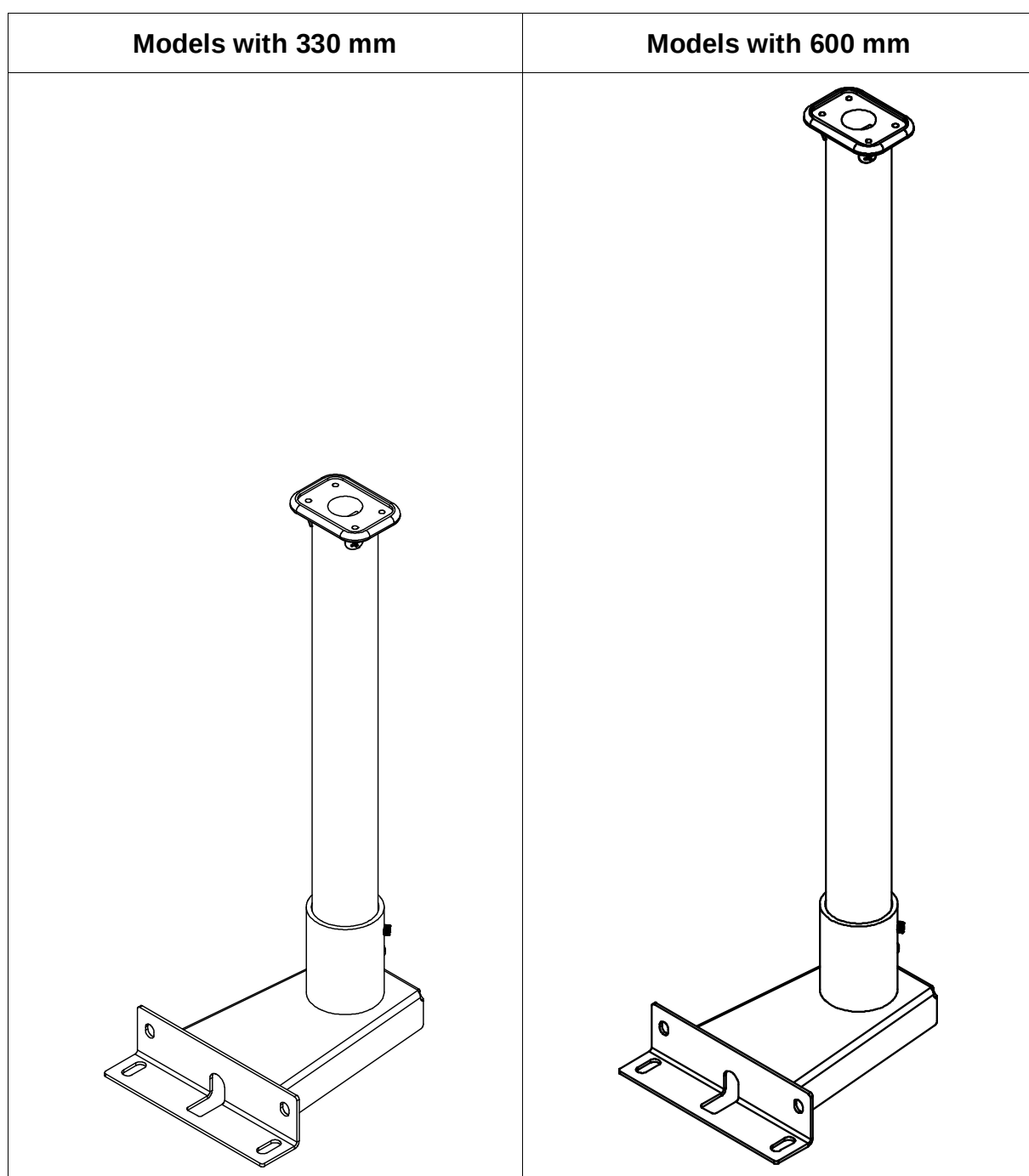









Models with 330 mm:

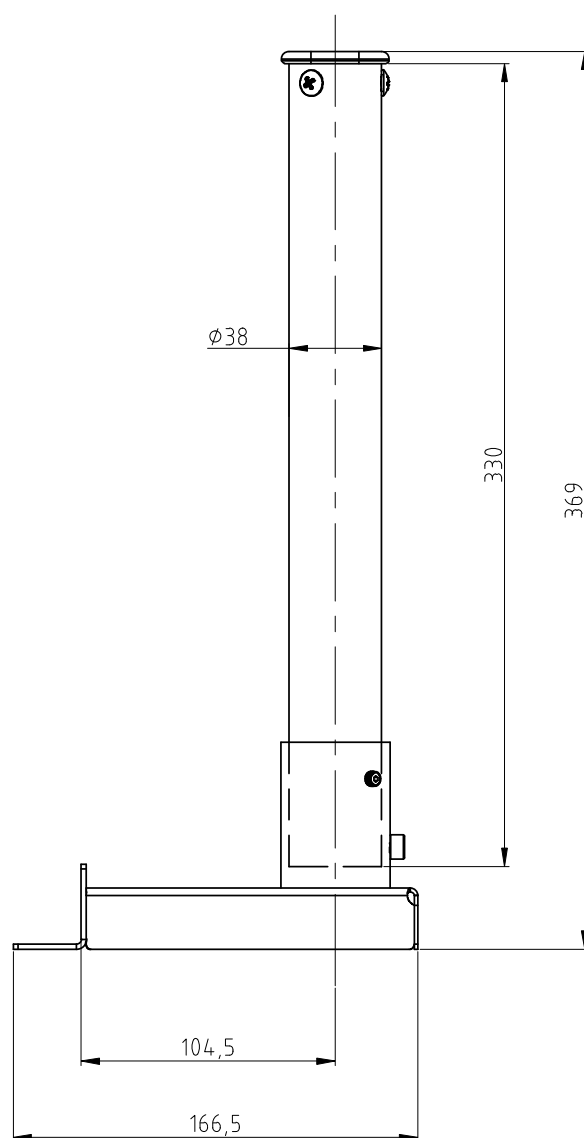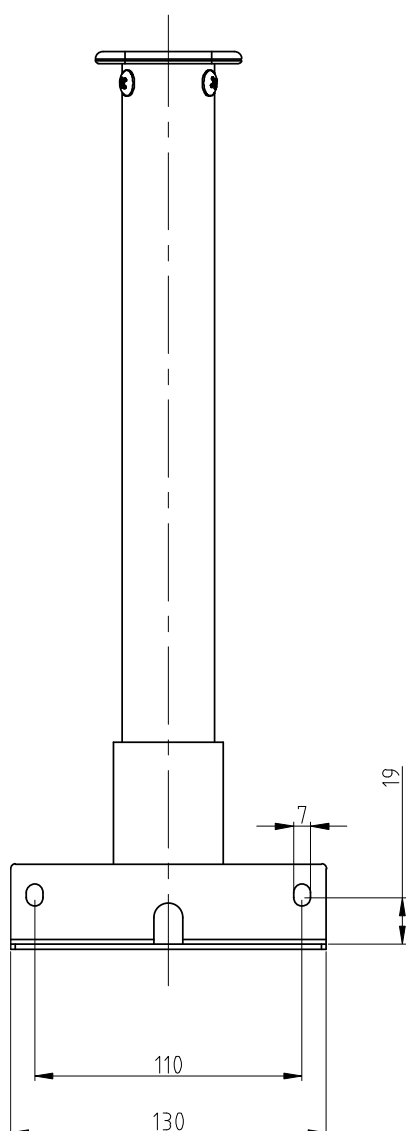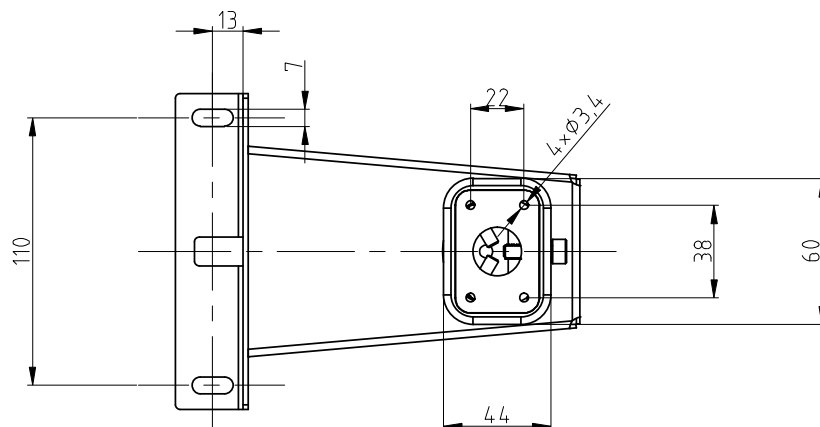

Models with 600 mm:

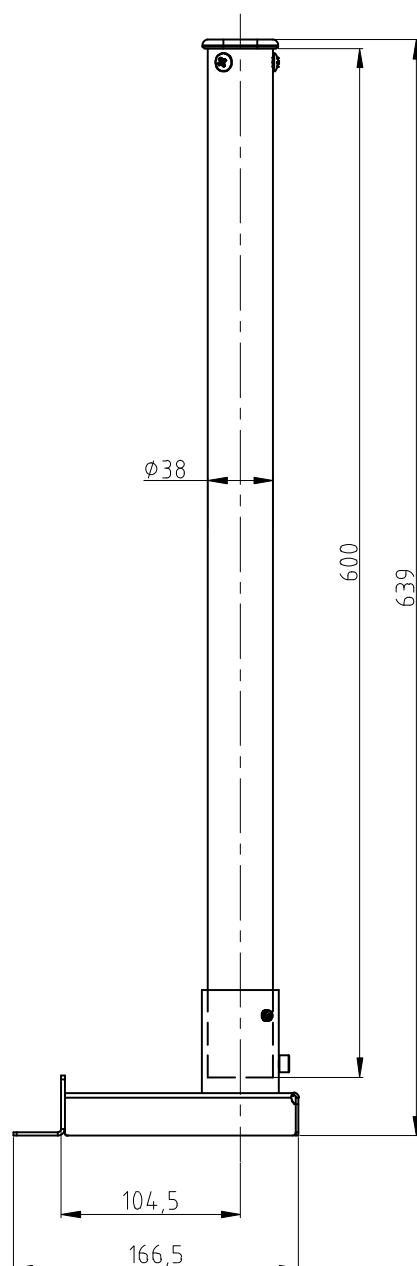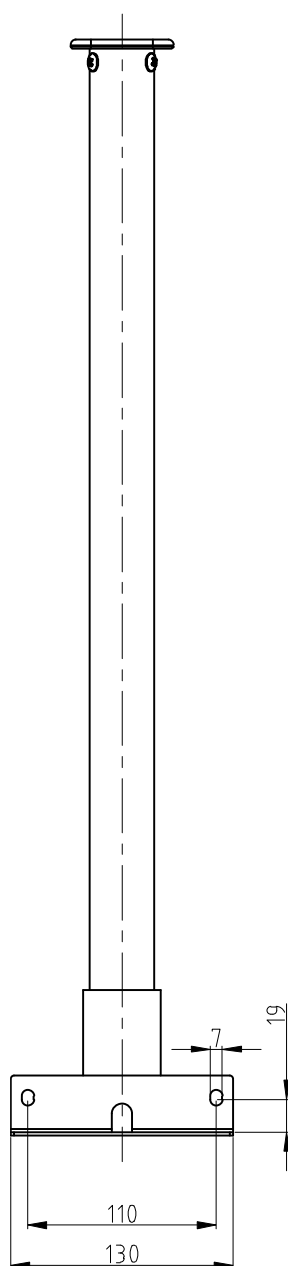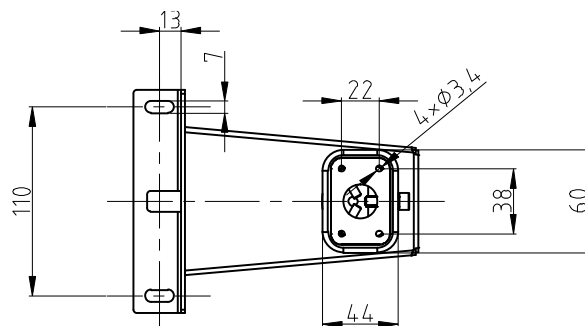

2 Scope of delivery

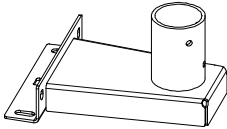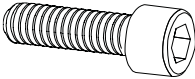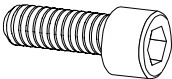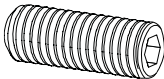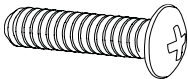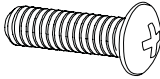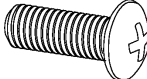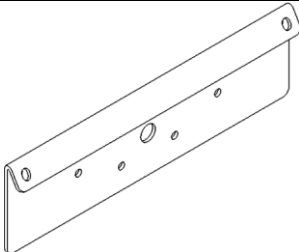
 A Tripod base 1 x	 B Tripod tube 1 x	 C Adapter 1 x
 S1 M6 x 16 4 x	 S2 Spring washer 4 x	 S3 Washer 4 x
 S4 M6 x 12 1 x	 S5 M6 x 10 2 x	 S6 M5 x 16 3 x
 S7 M4 x 10 4 x	 S8 M4 x 6 2 x	 S9 Adapter plate 1 x

3 Installation

Attach the stand base to the platform:


1. Position the stand base **(A)** centrally over the holes on the side of the platform.
2. First place the spring washer **(S2)** and then the washer **(S3)** on the screw **(S1)** (4 x).
3. Tighten the tripod base.

Make sure that you do not pinch or damage any cables.

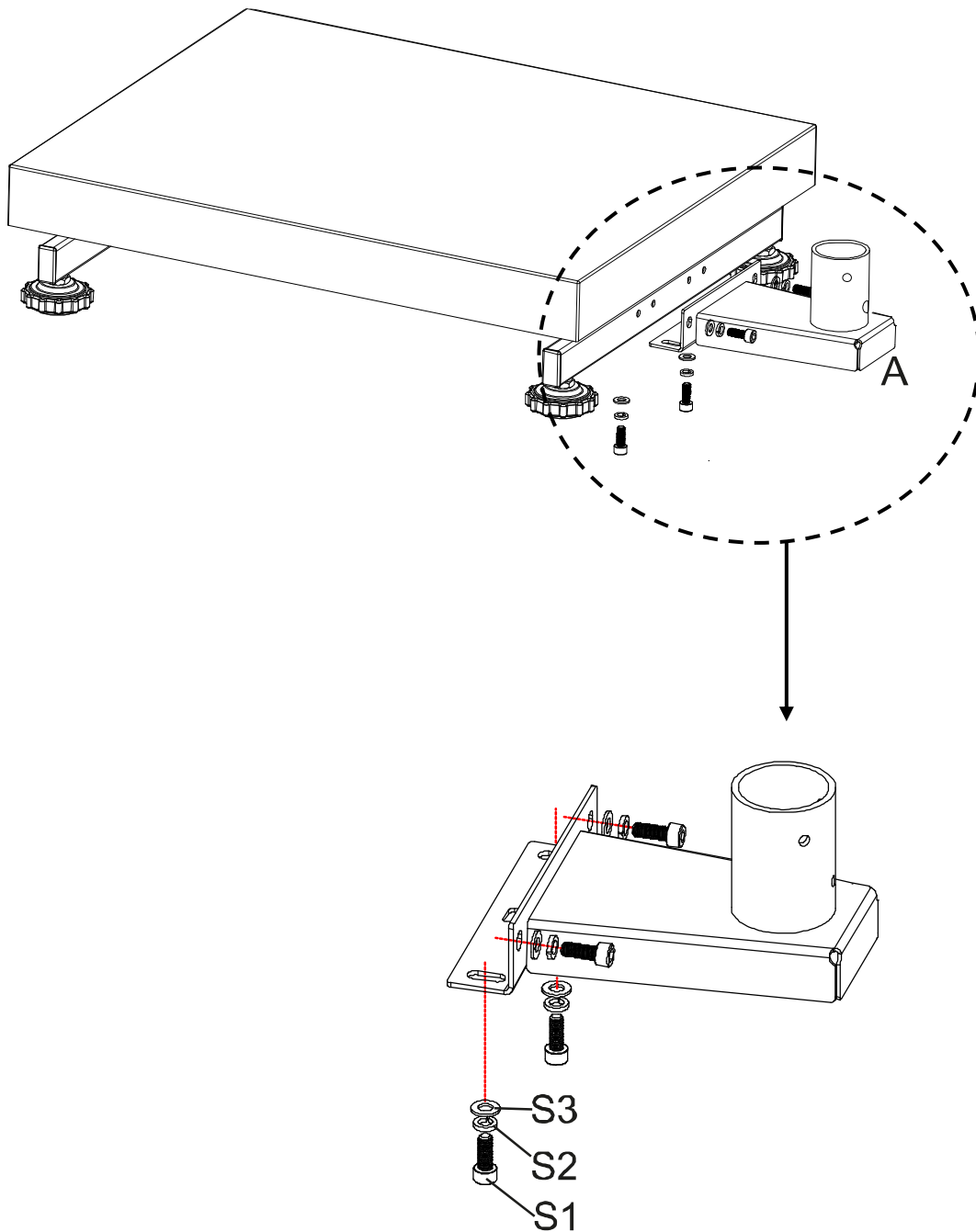


Attach the tripod tube to the tripod base:

4. Unscrew screws **S4** and **S5** from the stand tube.
5. Insert **the** tripod tube (**B**) into the tripod base. Ensure that the hole in the stand tube matches the hole **S4**.
6. Screw in screws **S5** again and fix the stand tube with screw **S4**.

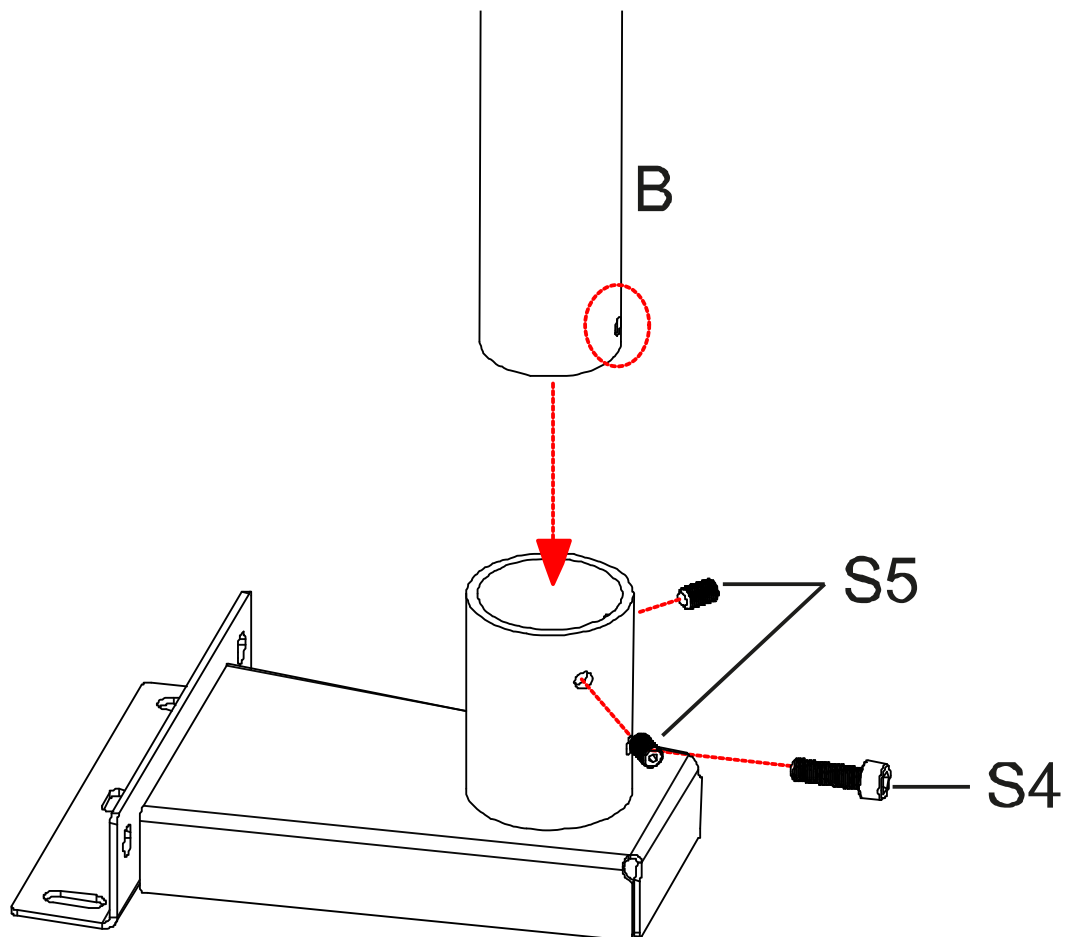

Attach the display unit to the stand tube:

7. Insert the adapter (**C**) into the stand tube (**B**) and fasten with the screws **S6**.
8. Attach the adapter plate **S9** to the holder for the display unit using the screws **S7**.

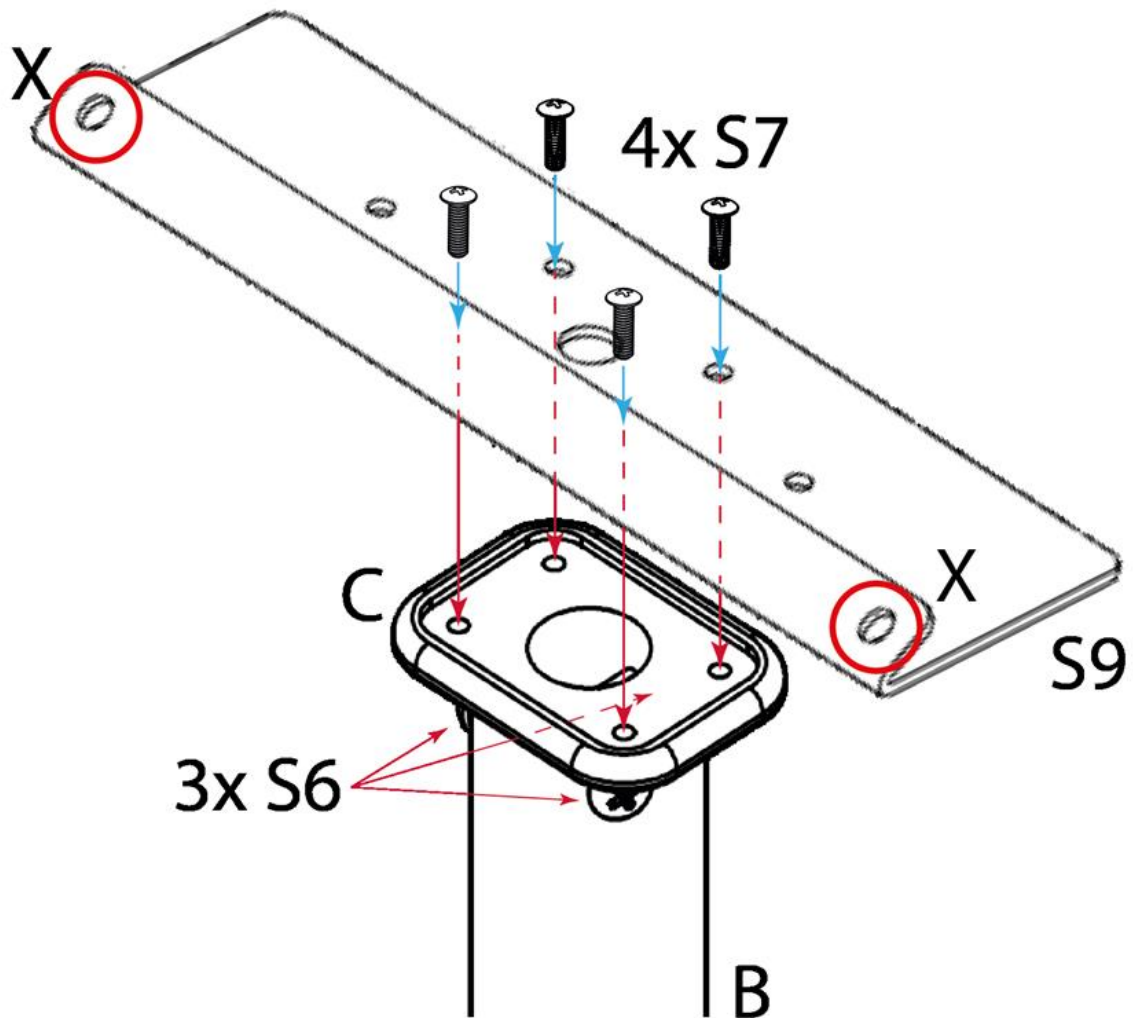

9. Fasten the display unit with the **S8** screws in the **X** holes.

4 Appendix

