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Ziegelei 1
D-72336 Balingen
E-Mail: info@kern-sohn.com

Tel: +49-[0]7433- 9933-0
Fax: +49-[0]7433-9933-149
Internet: www.kern-sohn.com

D

**Installationsanleitung
Akku**

Seite 2

GB

**Installation Information
Rechargeable battery**

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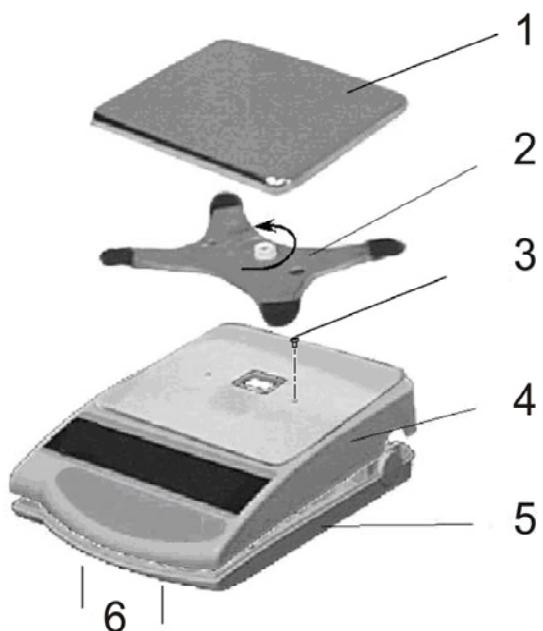
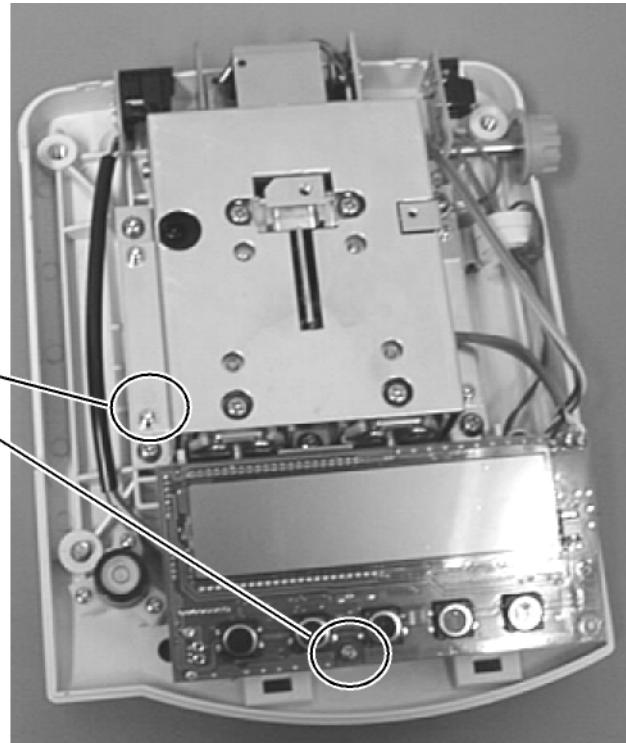
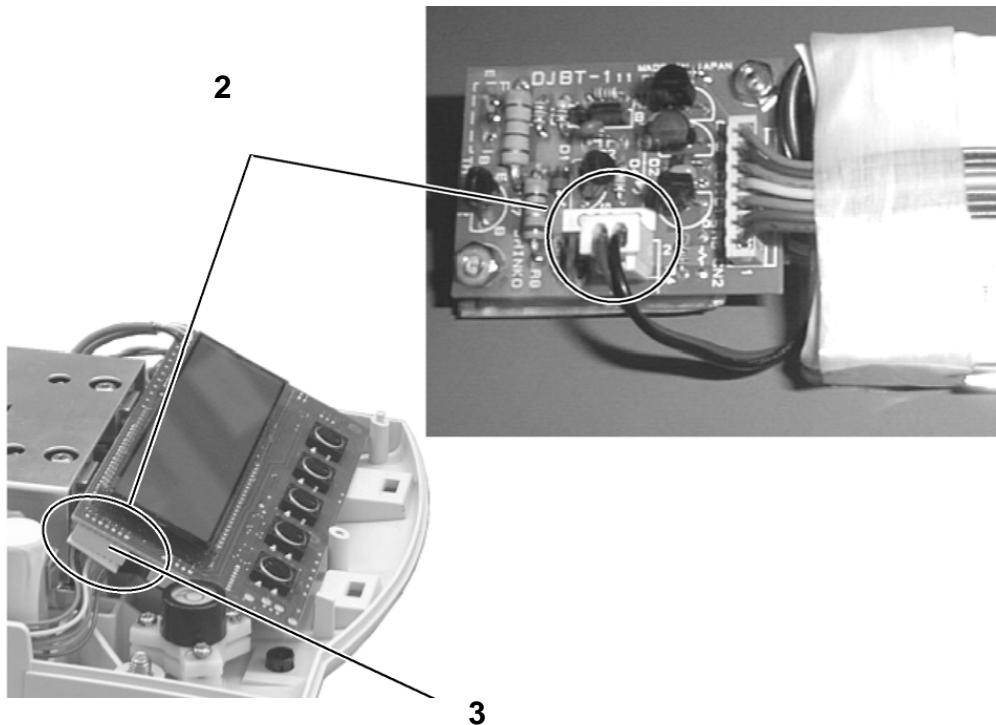
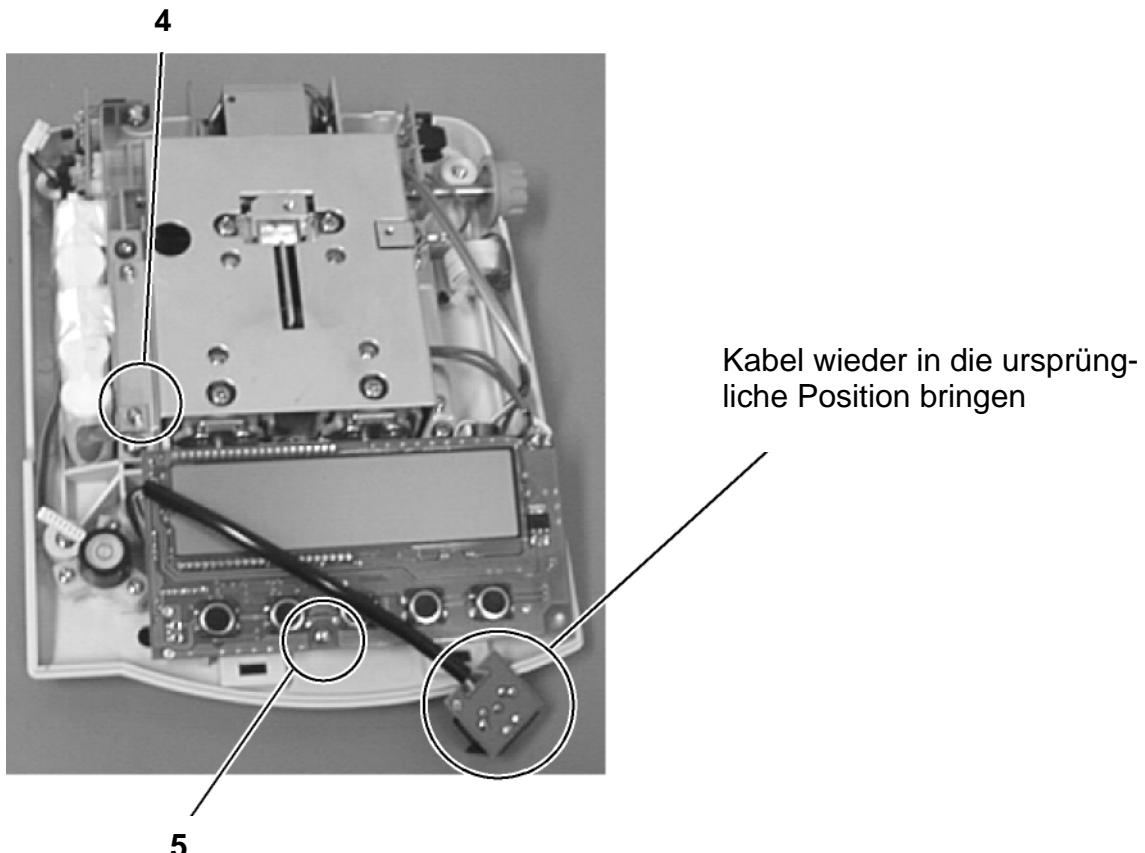
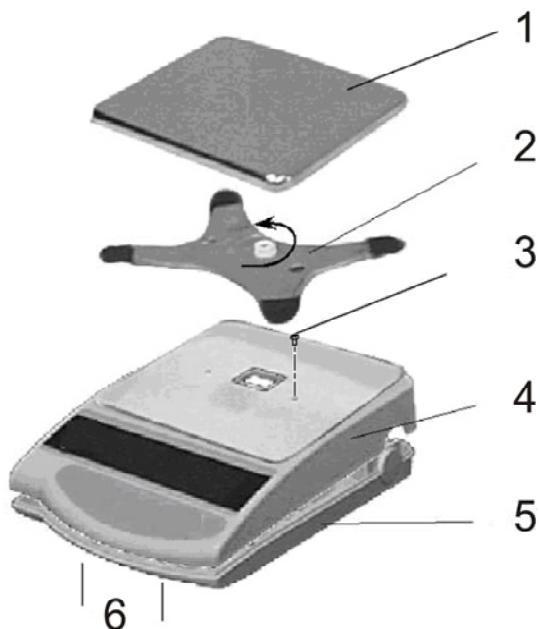
Installation Instructions

Rechargeable battery

KERN EG-A04 for models with max $\leq 820 \text{ g}$

How to remove the case:


- Disconnect the balance from the mains
- Remove weighing plate (1) and carrier of weighing plate (2)
- Remove both screws to the left and right of the bracket guide (3)


- Release both retaining hooks (6) on the lower housing section (5) and carefully remove the upper housing section (4) to the rear (please observe the upper housing section guides on the back of the balance).

To install rechargeable battery:

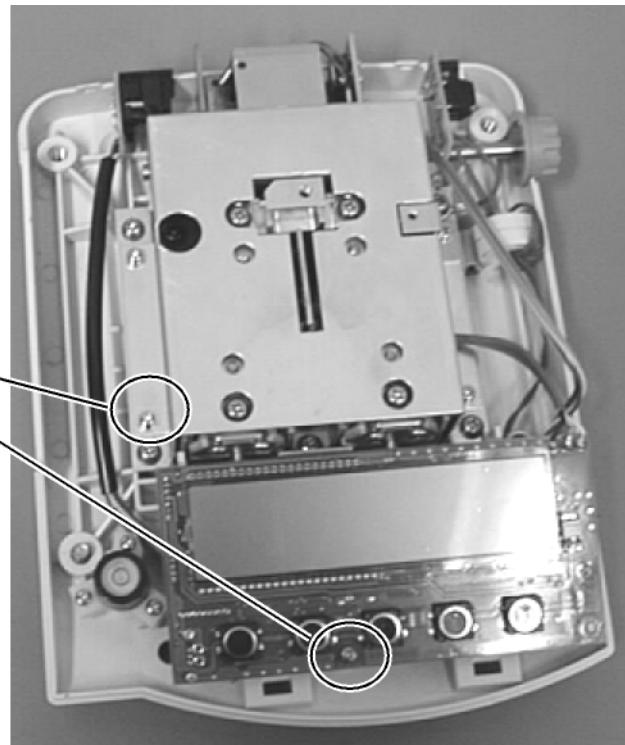
- Loosen and remove both fastening screws (1)

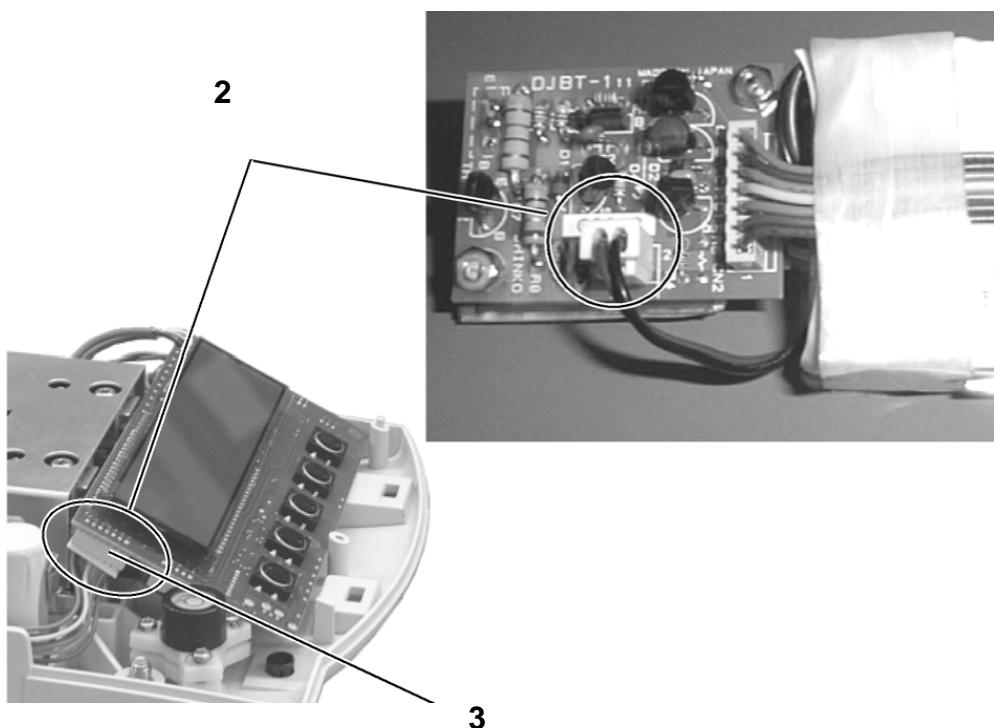- Remove the rechargeable battery from the packaging and begin by connecting the electric power supply to the circuit board of the rechargeable battery (2). Subsequently set up the plug-in connection to the computer circuit board (3) of the balance (CN5).

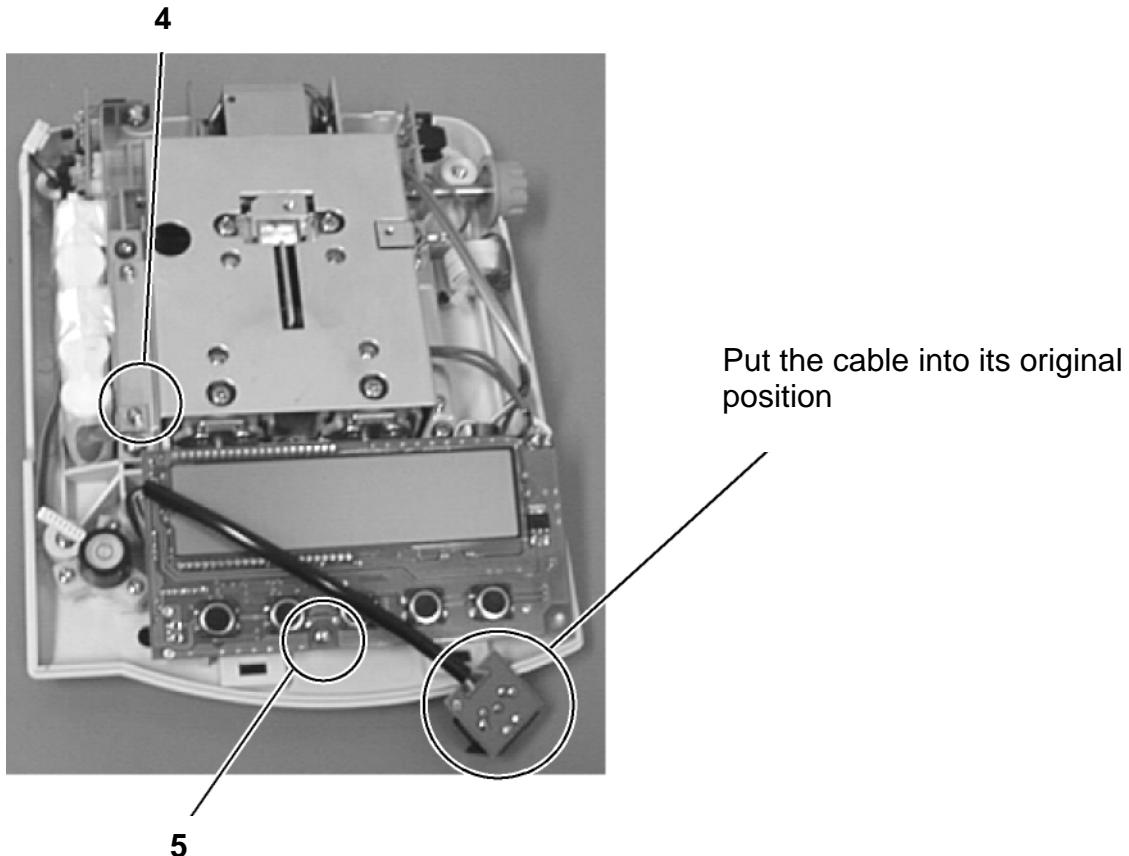- The rechargeable battery is placed on the left side of the housing in such a way that it can be screwed onto the balance through the existing mounting using the screw (4) loosened previously. Press lightly into the housing prior to this, (there is only one correct installation possibility). Now also re-attach the display using the loosened screw (5).

- Place the upper housing section on the rear guides and tilt forward until both retaining hooks on the lower housing section audibly lock into place. Screw down both screws to the left and right of the bracket guide tightly and re-secure the bracket. Attach the weighing plate.

Information:

Though the rechargeable battery is immediately operative, it should be charged for at least 8 hours using the mains adaptor before being used for the first time.

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