



KERN & Sohn GmbH

Ziegelei 1

D-72336 Balingen

E-Mail: info@kern-sohn.com

Phone: +49-[0]7433- 9933-0

Fax: +49-[0]7433-9933-149

Internet: www.kern-sohn.com

Installation Instructions

Signal lamp

KERN CFS-A03

Version 1.5

2025-06

en




CFS-A03-IA-e-2515

Configuration

The signal lamp **CFS-A03** can be adapted for different applications. For this proceed as follows:

To change the connections, loosen the 4 screws of the base plate and remove the base plate.

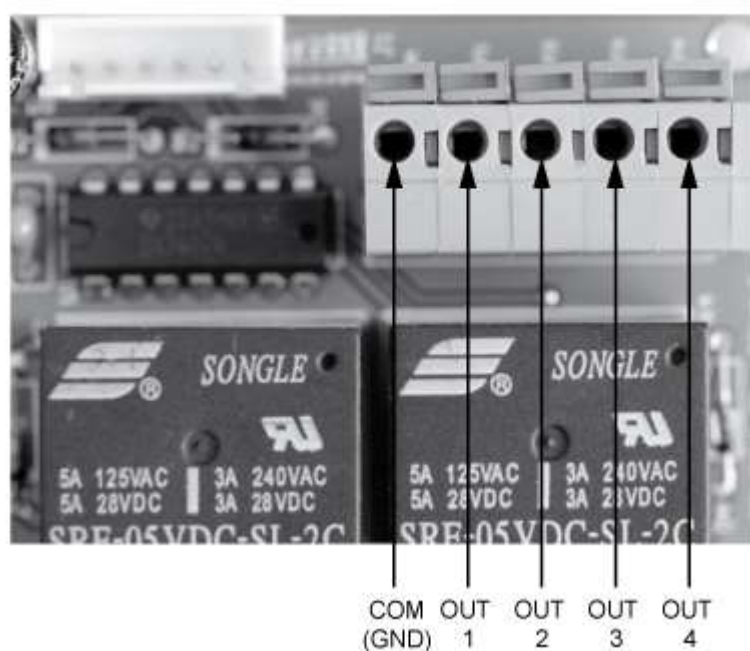

Do not alter connection „COM“ (GND)

Model GAB		Models KFB-TM, KFS-TM, CFS, CPB	
COM (GND)	Black	COM (GND)	Black
OUT 1	Yellow	OUT 1	Yellow
OUT 2	----	OUT 2	Red
OUT 3	Red	OUT 3	Green
OUT 4	Green	OUT 4	----
Models KUP-01 / KUM-01			
COM (GND)		Black	
OUT 1		Green	
OUT 2		Red	
OUT 3		Yellow	
OUT 4		----	