

Mounting the darkfield unit OBB-A1148 for the microscope series KERN OBE-1

The background to this is a technical modification of the condenser of the microscope, which causes a different use of the darkfield unit as well.

This circumstance is not yet included in the enclosed manual of the microscope.

Step 1:

Remove the bottom part of the condenser (see figure 1)



Figure 1

Step 2:

Screw the dark field unit supplied (see figure 2) into the lower part of the condenser as far as it will go.



Figure 2

Note:

To switch back to the brightfield method, you must attach the standard bottom part of the condenser anew.