



**KERN & Sohn GmbH**

Ziegelei 1  
72336 Balingen-Frommern  
Germany

+0049-[0]7433-9933-0  
+0049-[0]7433-9933-149  
info@kern-sohn.com

# Installation Instructions

## Interface Adapter with Cable for Extension Box

### KERN KUP-13

Type TYKUP-13-A  
Version 1.2  
2025-05  
GB



You will find the current version of these instructions also online under:  
<https://www.kern-sohn.com/shop/de/DOWNLOADS/>  
Under the column Operating instructions

TYKUP-13-A-IA-e-2512



**Interface Adapter with Cable for Extension Box**  
Version 1.2 2025-05  
**Installation Instructions**

---

---

## Contents

<b>1</b>	<b>Scope of delivery .....</b>	<b>2</b>
<b>2</b>	<b>General information.....</b>	<b>2</b>
<b>3</b>	<b>Installation .....</b>	<b>2</b>

### 1 Scope of delivery

- Interface Adapter with Cable for Extension Box
- Installation Instructions

### 2 General information

**Cable length:** 0,15 m

<b>i</b>	<ul style="list-style-type: none"><li>• Via the KERN Extension Box up to 3 interfaces of your choice can be connected. These can be used in parallel and simultaneously.</li><li>• Cascading is not possible!</li><li>• <b>Only KERN KUP adapters may be connected to the 15-pol-sub-D-plug of the device!</b></li></ul>
----------	--

### 3 Installation

1. Switch off the appliance
2. Plug in the interface adapter at the KUP connection
3. Switch on the appliance. The interface adapter is automatically recognized by the appliance. However, it may take up to one minute before the interfaces can be used.