



KERN & Sohn GmbH

Brickyard 1
D-72336 Balingen
E-mail: info@kern-sohn.com

Tel: +49-[0]7433- 9933-0
Fax: +49-[0]7433-9933-149
Internet: www.kern-sohn.com

Installation instructions Tripod

KERN IFB-A01 / A02

Version 1.3
2025-06
en



IFB-A01-A02-IA-e-2513



Tripod

Version 1.3 2025-06

Installation instructions

Table of contents




1	Technical data.....	3
1.1	Dimensions.....	4
2	Scope of delivery	6
3	Installation	7

1 Technical data




KERN	IFB-A01	IFB-A02
Tripod height (mm)	330	600
For platform sizes (mm)	300 x 240 400 x 300 500 x 400 650 x 500	400 x 300 500 x 400 650 x 500

1.1 Dimensions






Models with 330 mm:

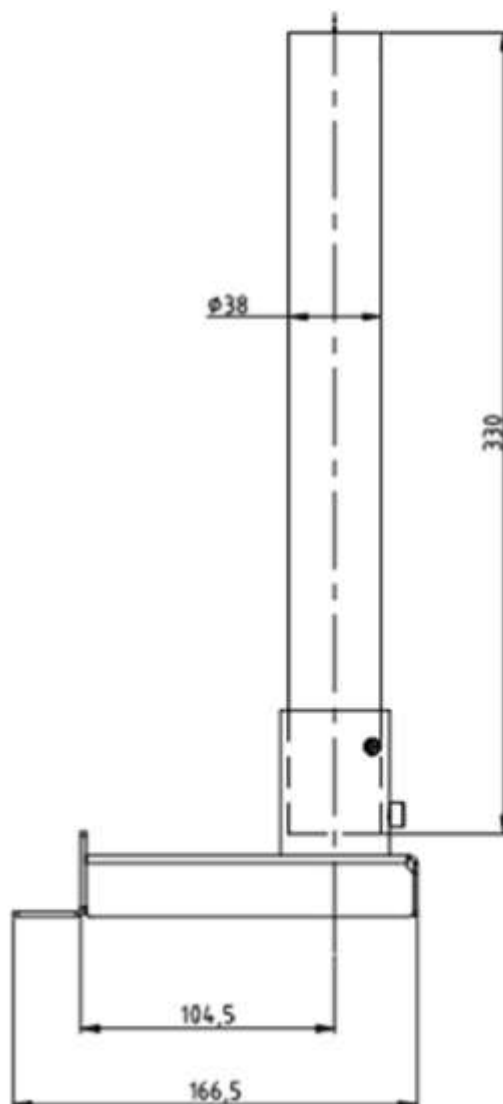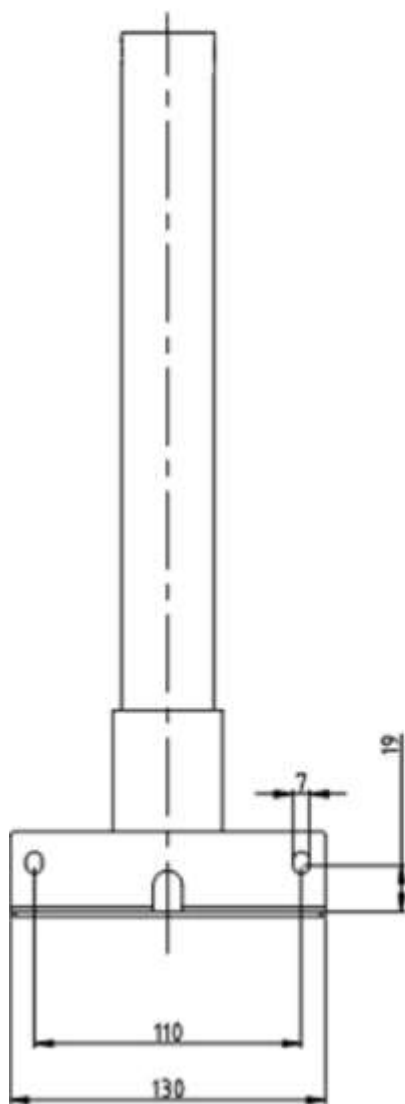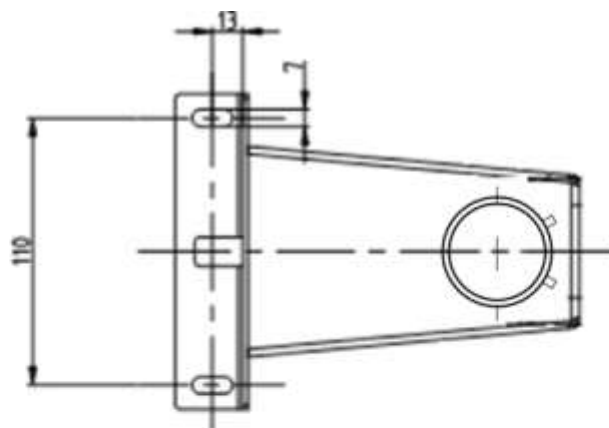

Models with 600 mm:

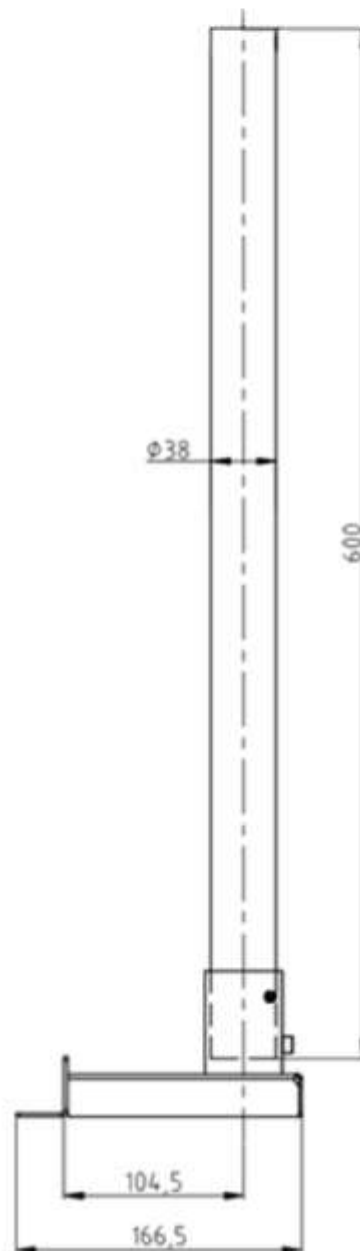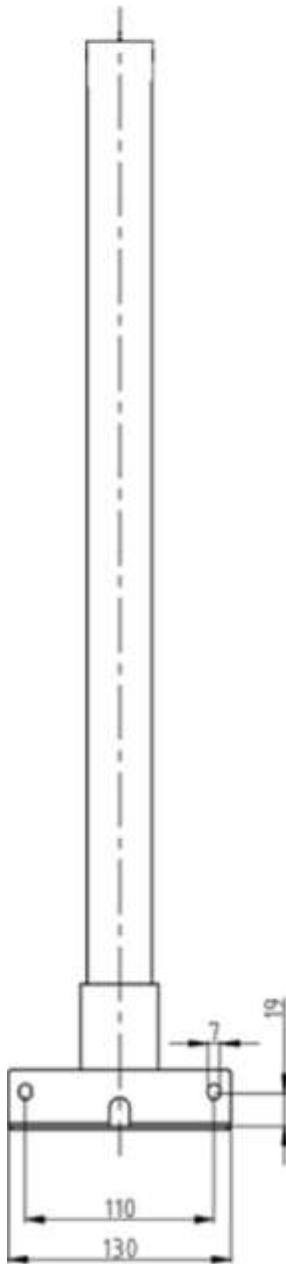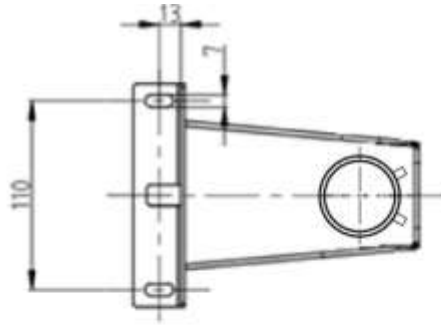

2 Scope of delivery

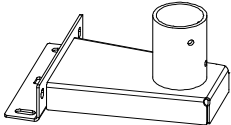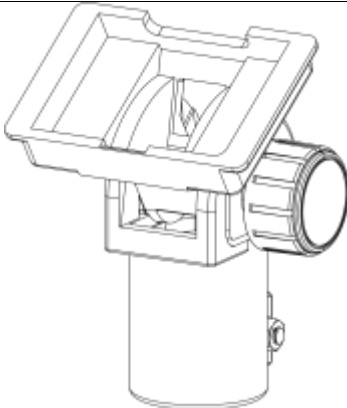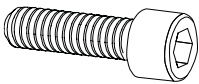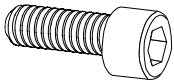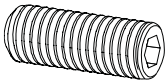 A Tripod base 1 x	 B Tripod tube 1 x	 C Holder for display unit 1 x
 S1 M6 x 16 4 x	 S2 Spring washer 4 x	 S3 Washer 5 x
 S4 M6 x 12 1 x	 S5 M6 x 10 2 x	 S6 M8 x 20 1 x
 S7 M8 nut 1 x		


C, **S3** (1x), **S6** and **S7** are supplied pre-assembled.

3 Installation

Attach the stand base to the platform:


1. Position the stand base **(A)** centrally over the holes on the side of the platform.
2. First place the spring washer **(S2)** and then the washer **(S3)** on the screw **(S1)** (4 x).
3. Tighten the tripod base.

Make sure that you do not pinch or damage any cables.

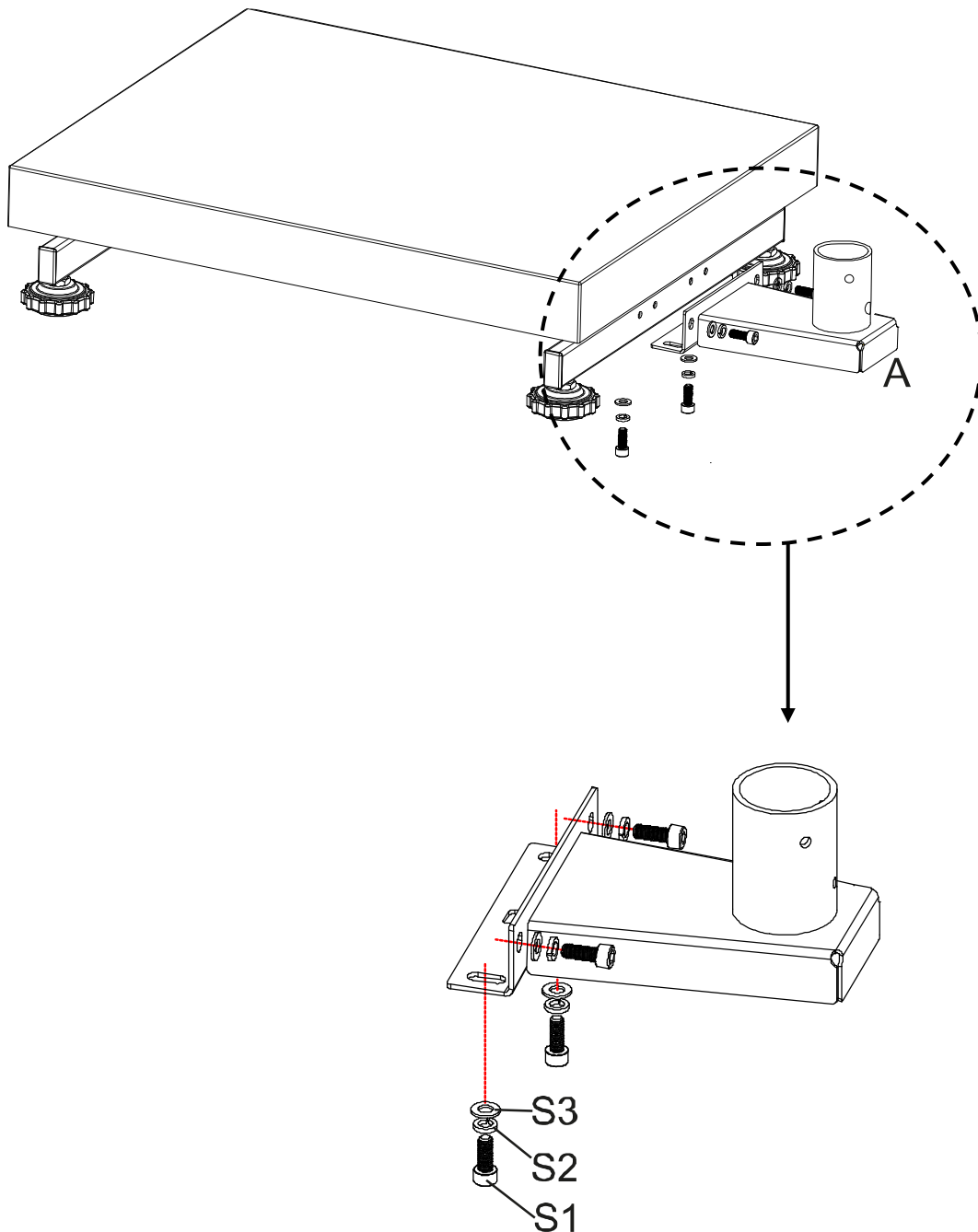


Attach the tripod tube to the tripod base:

4. Unscrew screws **S4** and **S5** from the stand tube.
5. Insert **the** tripod tube (**B**) into the tripod base. Ensure that the hole in the stand tube matches the hole **S4**.
6. Screw in screws **S5** again and fix the stand tube with screw **S4**.

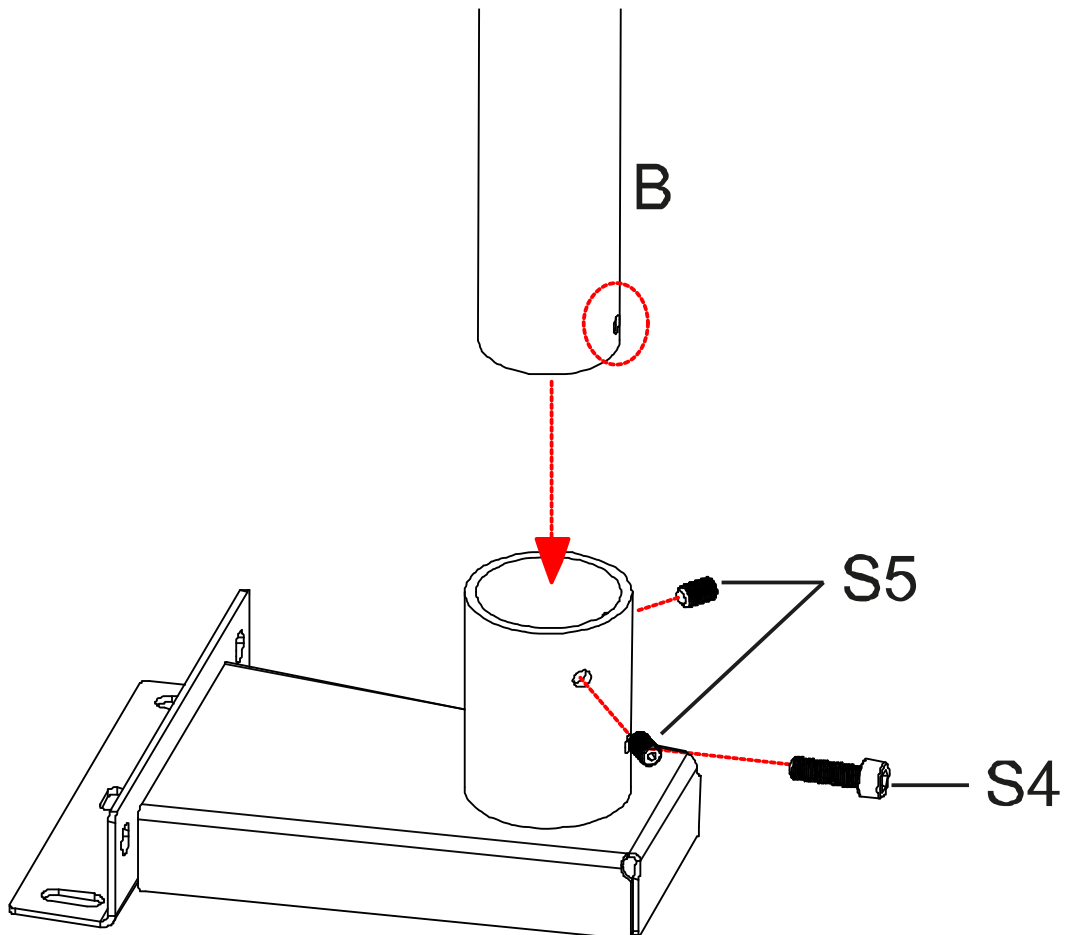

Attach the display unit to the stand tube:

7. Insert the holder for the display unit (**C**) into the stand tube and fasten with the screws **S6**.

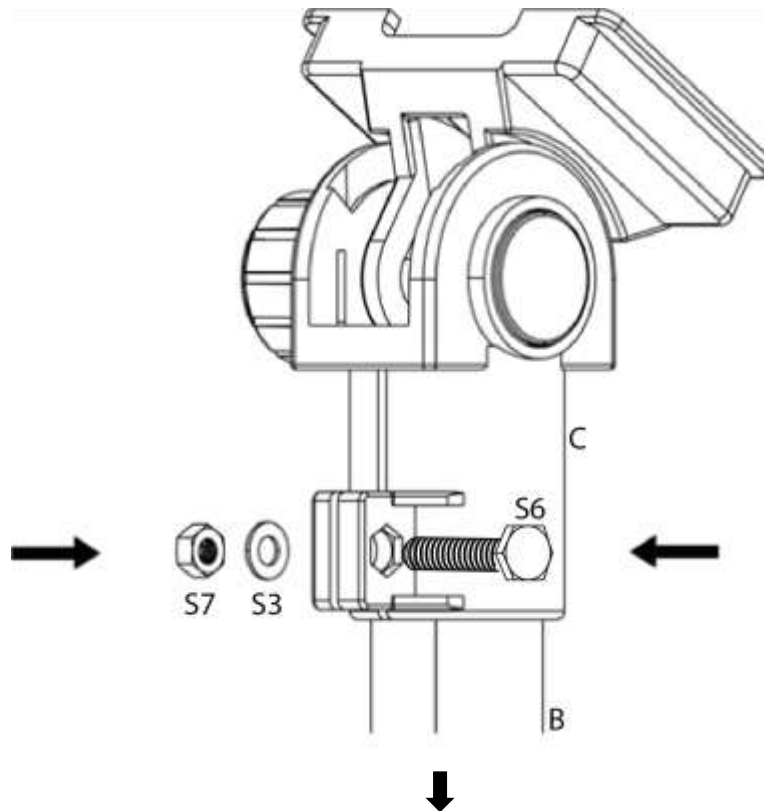

⇒ Insert the holder (**C**) for the display unit into the upper end of the stand tube (**B**) and secure with the screw (**S6**).

Counter on the opposite side with the washer (**S3**) and the nut (**S7**).