



KERN & Sohn GmbH

Ziegelei 1
D-72336 Balingen
E-Mail: info@kern-sohn.com

Tel: +49-[0]7433-9933-0
Fax: +49-[0]7433-9933-149
Internet: www.kern-sohn.com

Operating Instructions KERN EasyTouch

EasyTouch Installation setup User manual

Version 1.1
2022-09
GB



Contents


1.0 Set up of KERN Easy Touch	3
1.1 Download and Installation	3
1.2 Kern account activation	4
2.0 User without a registered Kern account	4
2.1 Register a KERN account	4
2.2 Enable the trial activation	7
2.2.1 Trial activation	7
2.3 Admin creation	10
2.4 Purchase licenses	13
2.5 License activation	14
3.0 User with the registered Kern account	19

1.0 Set up of KERN Easy Touch


A proper user account is required to download the application from any store. The Basic ET – License [SET – 01] is always required since it’s the base for working with KERN Easy-Touch.

1.1 Download and Installation


- Go to the Microsoft store and search for the Kern Easy Touch 2.0 in the search bar.

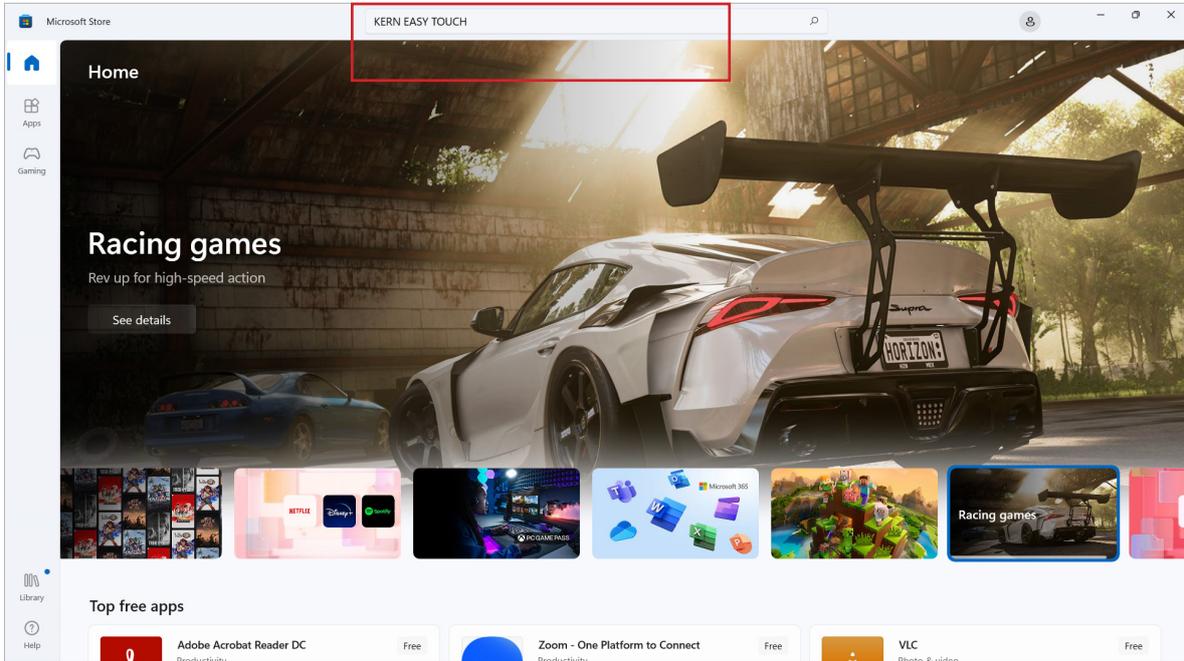

- Then there you can Easy Touch 2.0 application which is ready to install. There click on “install”

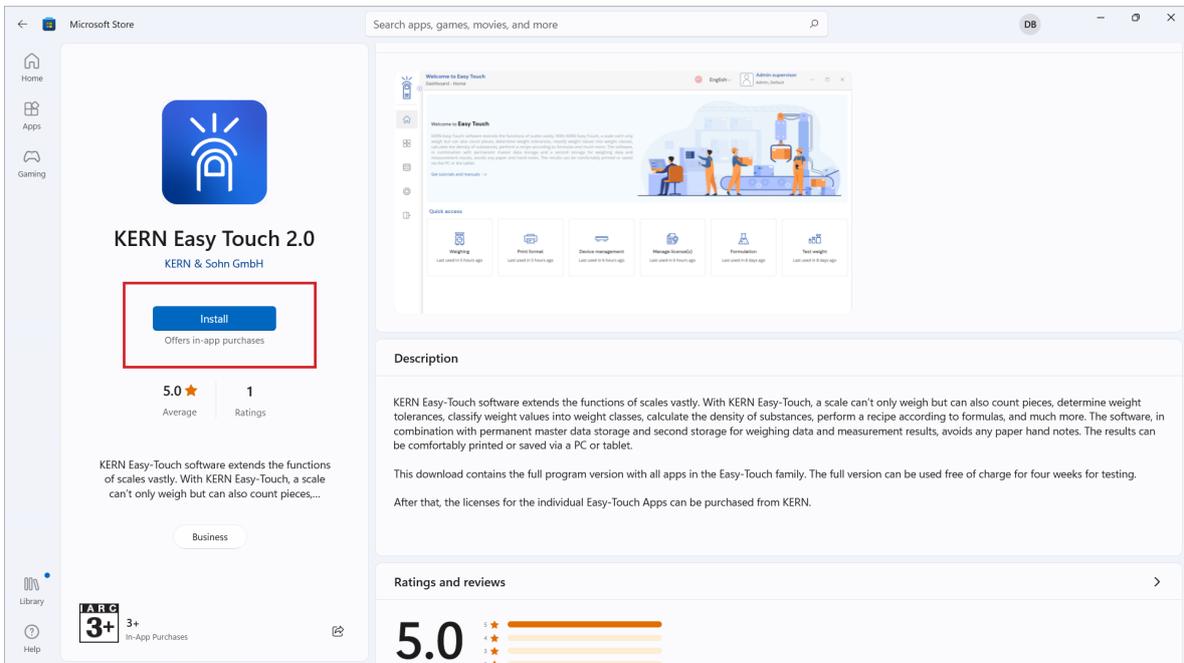

- When you click on install, the application will be downloaded and ready for use. Clicking on open will launch the application.

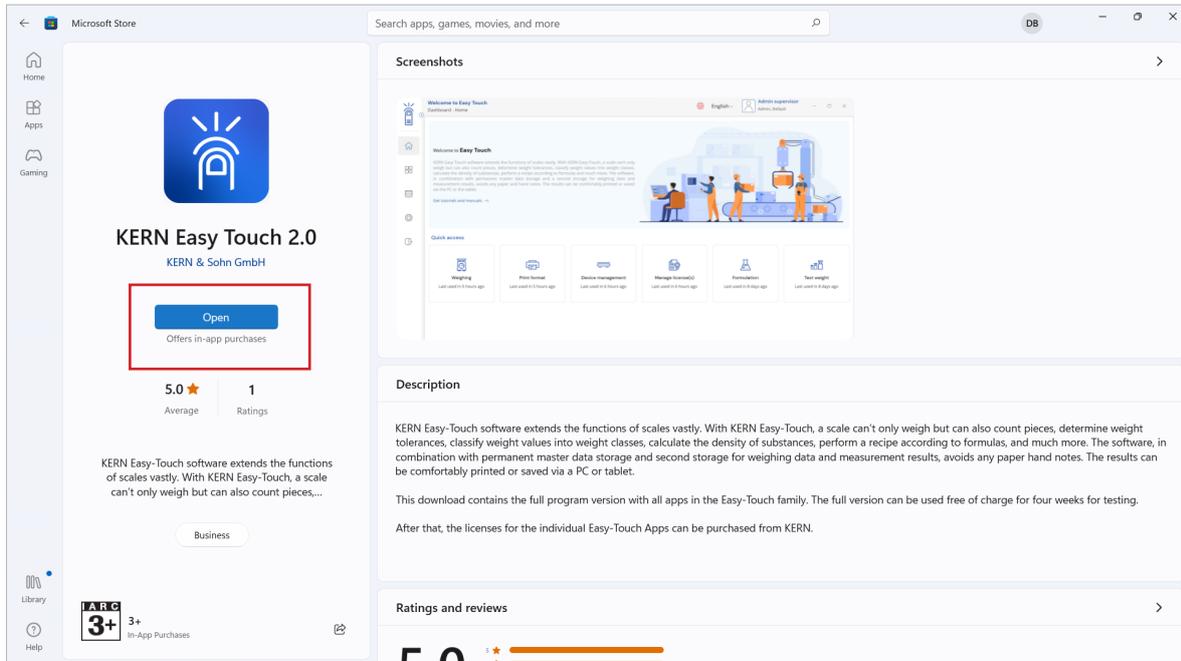

1.2 Kern account activation

KERN account registration is mandatory to utilise the license of Easy Touch.


2.0 User without a registered Kern account

The following steps has to be performed for using the Easy Touch functions for the user without a KERN account.


- [Register a KERN account](#)
- [Enable the trial license \(or purchase\)](#)
- [Admin creation for your system](#)
- [Purchase the functions](#)
- [Activation of purchased licenses](#)

2.1 Register a KERN account


The user can click the register KERN account hyperlink at the bottom left of the “Login and activate” screen to register after opening the Easy Touch application.

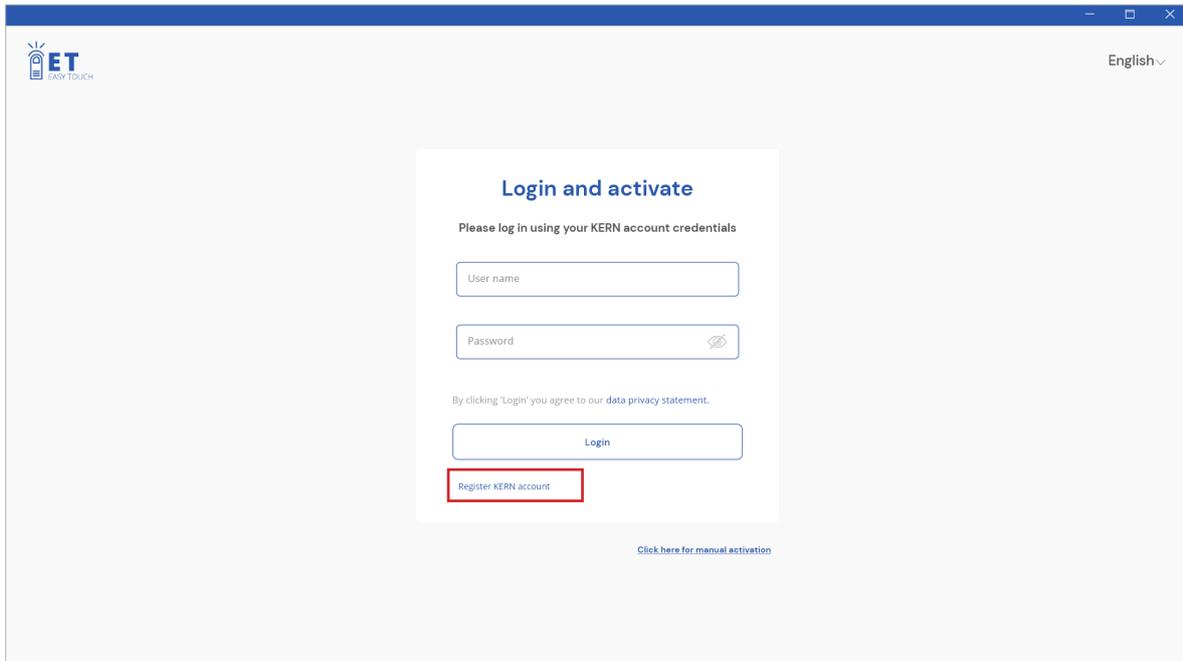

- After clicking the create account link the user will be navigated to the kern registration screen.

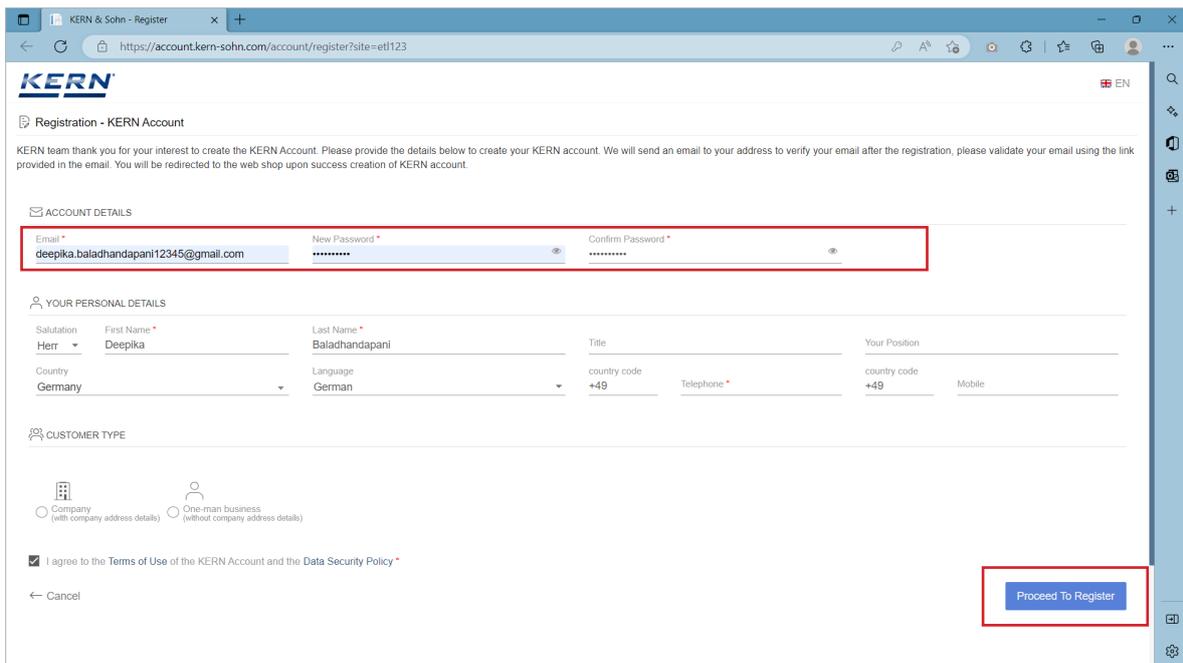

- Fill out and the mandatory fields and click the “proceed to register” button at the bottom of the screen.
- After proceeding to register, you will receive a confirmation message.

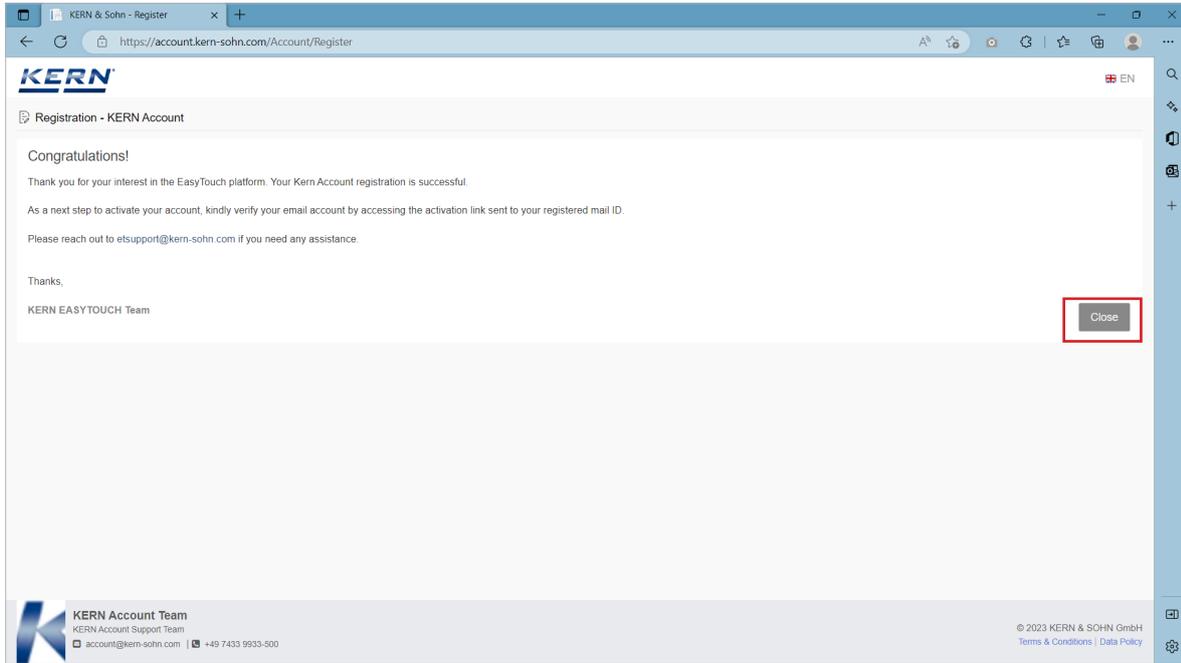


- An auto generated email will be sent to the registered email address with an activation link enclosed.
- Click on the hyperlink for verification. As soon as the email address is verified, the registration has been finished successfully.




- Now, login the Easy Touch application with the registered mail id and password

English

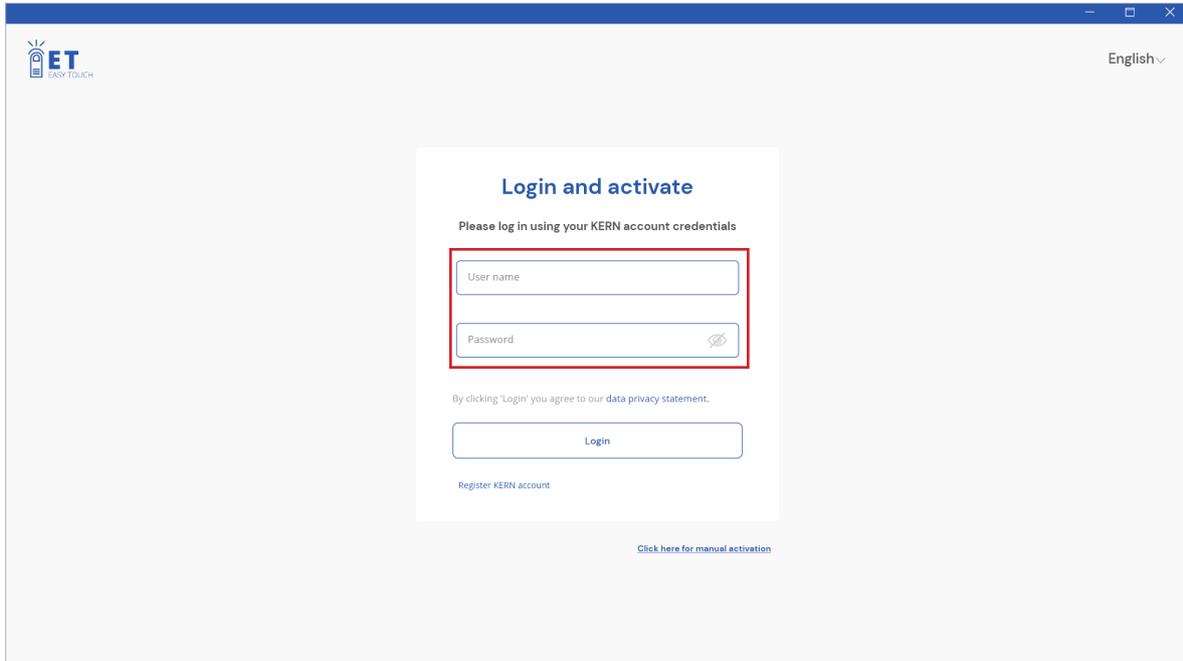

2.2 Enable the trial activation

- Activate the trial version by clicking on “start trial”.

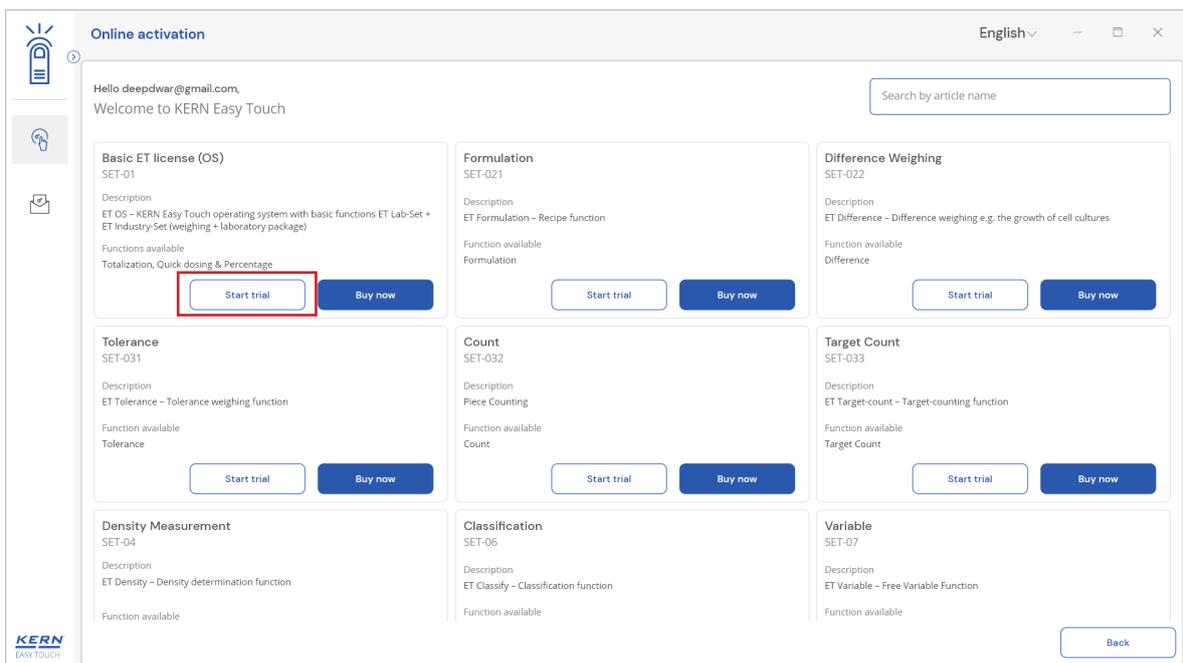

- Kindly note, the trial for the particular article will be applicable for the mentioned days from the date of activation and it will automatically deactivated once the trial is over.

2.2.1 Trial activation

- Click on the required license to start the trial activation.

Kindly note, the activation of basic ET license (SET 01) is mandatory for other article activation.

Online activation

Hello deepdwar@gmail.com,
Welcome to KERN Easy Touch

Search by article name

Basic ET license (OS) SET-01 Description ET OS - KERN Easy Touch operating system with basic functions ET Lab-Set + ET Industry-Set (weighing + laboratory package) Functions available Totalization, Quick dosing & Percentage Start trial Buy now	Formulation SET-021 Description ET Formulation - Recipe function Function available Formulation Start trial Buy now	Difference Weighing SET-022 Description ET Difference - Difference weighing e.g. the growth of cell cultures Function available Difference Start trial Buy now
Tolerance SET-031 Description ET Tolerance - Tolerance weighing function Function available Tolerance Start trial Buy now	Count SET-032 Description Piece Counting Function available Count Start trial Buy now	Target Count SET-033 Description ET Target-count - Target-counting function Function available Target Count Start trial Buy now
Density Measurement SET-04 Description ET Density - Density determination function Function available	Classification SET-06 Description ET Classify - Classification function Function available	Variable SET-07 Description ET Variable - Free Variable Function Function available

Basic set (SET-01) is needs to be purchased/activated

- Click on the start trial for the required functions

Online activation


Hello deepdwar@gmail.com,
Welcome to KERN Easy Touch

Search by article name


Basic ET license (OS) SET-01 Description ET OS - KERN Easy Touch operating system with basic functions ET Lab-Set + ET Industry-Set (weighing + laboratory package) Functions available Totalization, Quick dosing & Percentage Start trial Buy now	Formulation SET-021 Description Formulation Function available Formulation Start trial Buy now	Difference Weighing SET-022 Description Difference Weighing Function available Difference Start trial Buy now
Tolerance SET-031 Description Check weighing Function available Tolerance Start trial Buy now	Count SET-032 Description Piece Counting Function available Count Start trial Buy now	Target Count SET-033 Description Target Count Function available Target Count Start trial Buy now
Density Measurement SET-04 Description Density Measurement Function available	Classification SET-06 Description Classification Function available	Variable SET-07 Description Variable Function available

Back


English

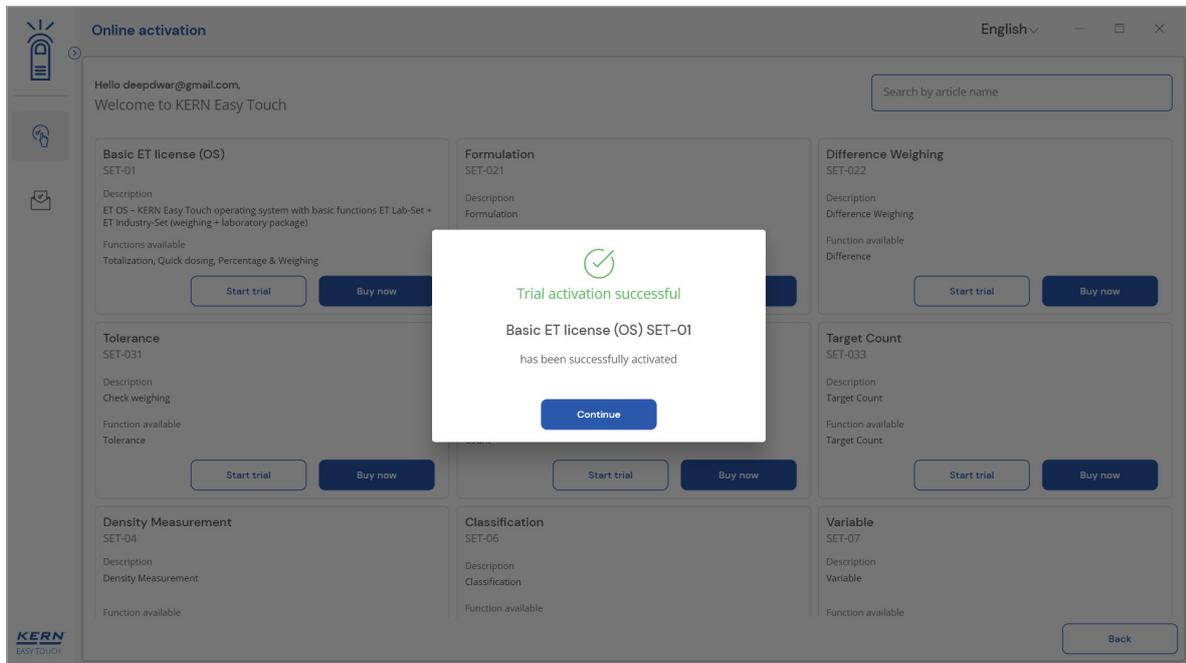

- The trial period would be mentioned according to the function and user will be allowed to purchase the function.

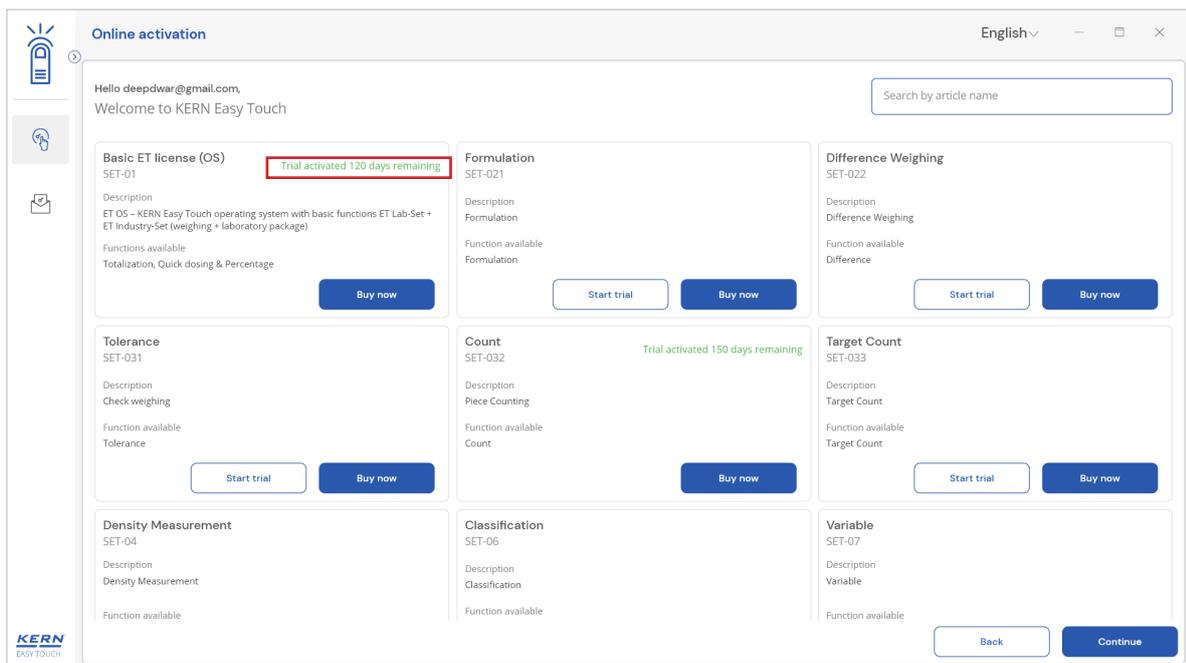


- User can activate the trial for required functions and click on continue to proceed in using the functions

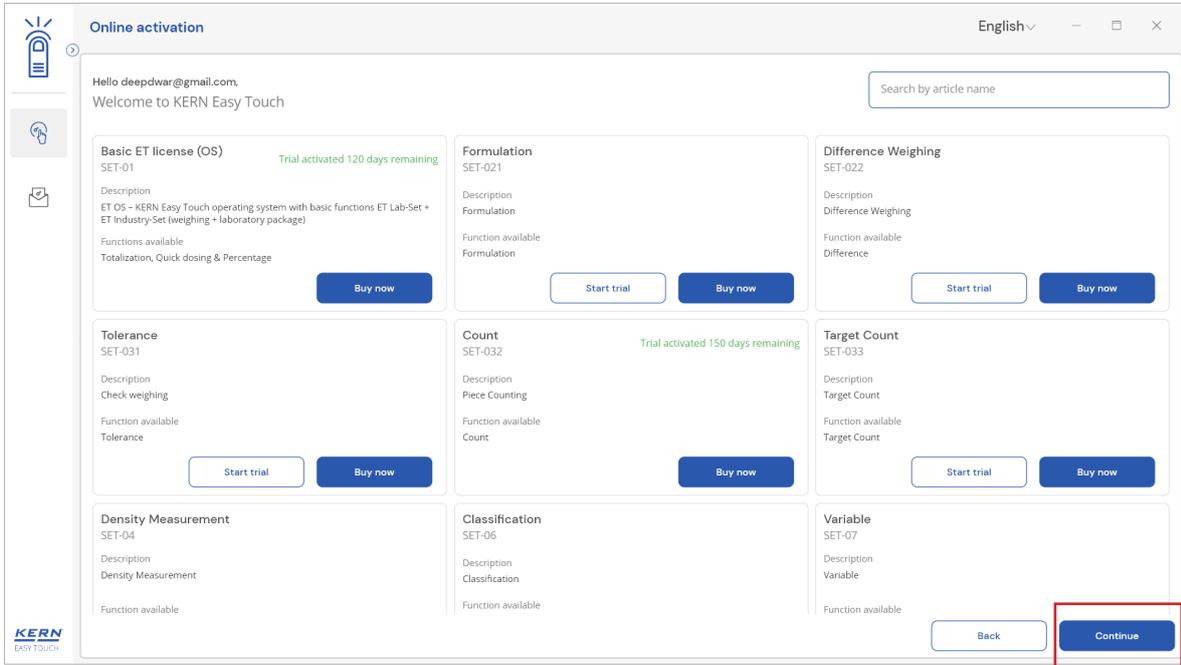

- Create your admin account and login and you are set up to utilise the activated functions.

2.3 Admin creation

When you log in to the Easy Touch application for the first time, you will be prompted to create an admin user for the Easy Touch application post the activation of licenses.

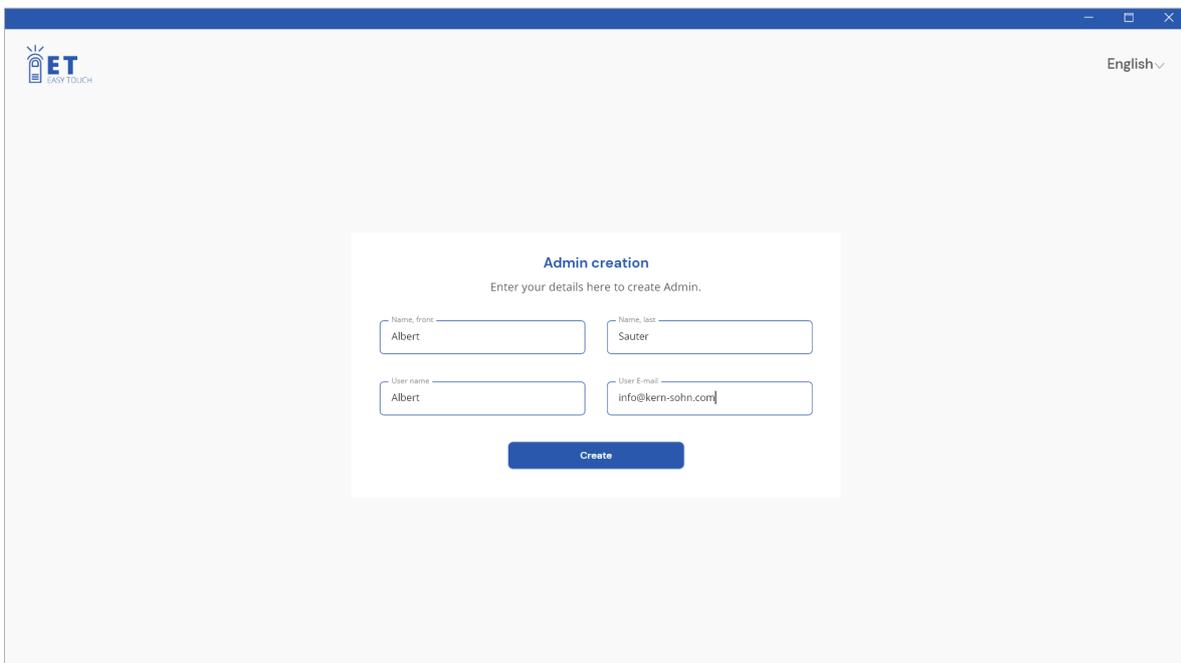

- Fill all the required details and click the create

ET EASY TOUCH English

Admin creation
Enter your details here to create Admin.

Name, first: Albert
Name, last: Sauter
User name: Albert
User E-mail: info@kern-sohn.com

Create

- A temporary password will be generated after clicking the create button and it will be displayed. Additionally, you will receive this temporary password via email to the mail ID that you have specified in the admin creation screen.


ET EASY TOUCH English

Admin creation
Please use the temporary password to log in

Temporary password: 945207

Copy & proceed

- Now upon clicking on “copy and proceed” the otp will be copied and then the user login screen appears.
- Now enter the created username and the temporary password and click on login.

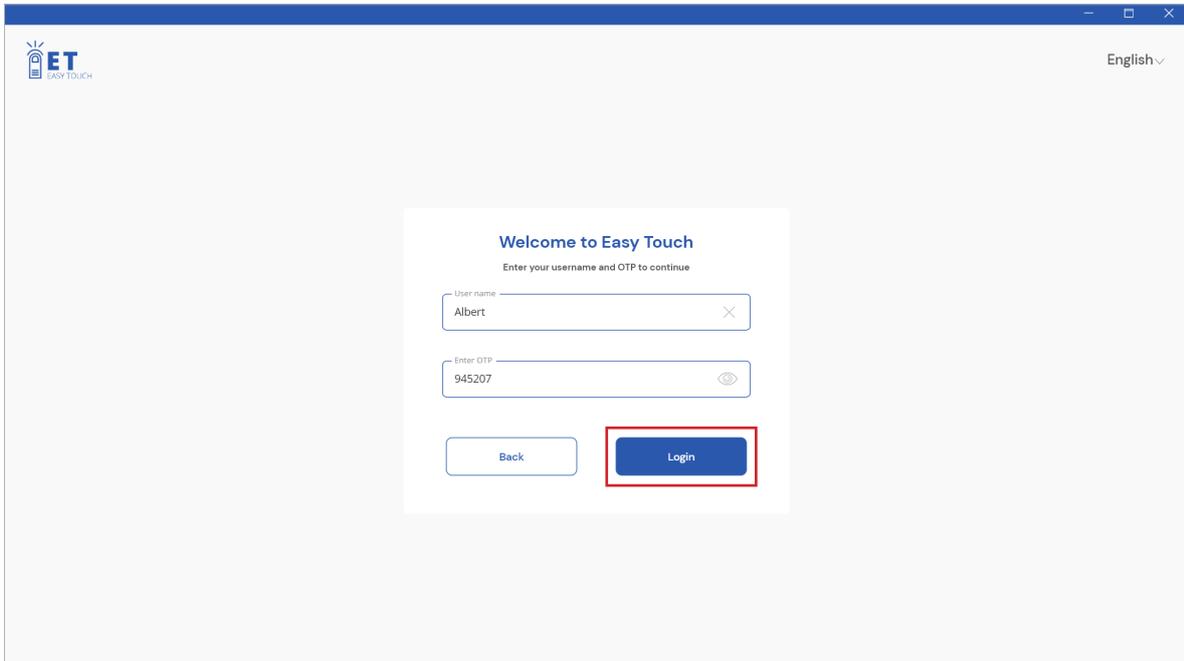

ET
EASY TOUCH

English

Welcome to Easy Touch


Enter your username and OTP to continue

User name
Albert

Enter OTP
945207

Back Login

- Reset the password and click on next. Your admin account is now created and new password is reset.

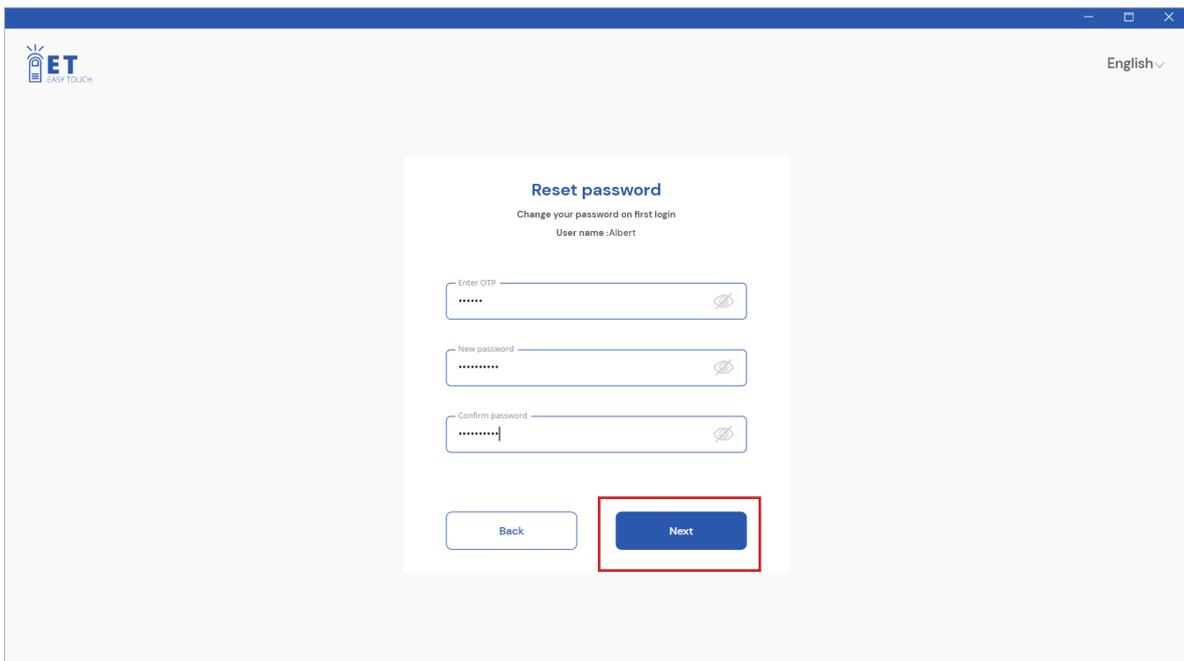

ET
EASY TOUCH

English

Reset password

Change your password on first login
User name :Albert


Enter OTP
.....

New password
.....

Confirm password
.....

Back Next

- Enter the user name and new password and click on login.

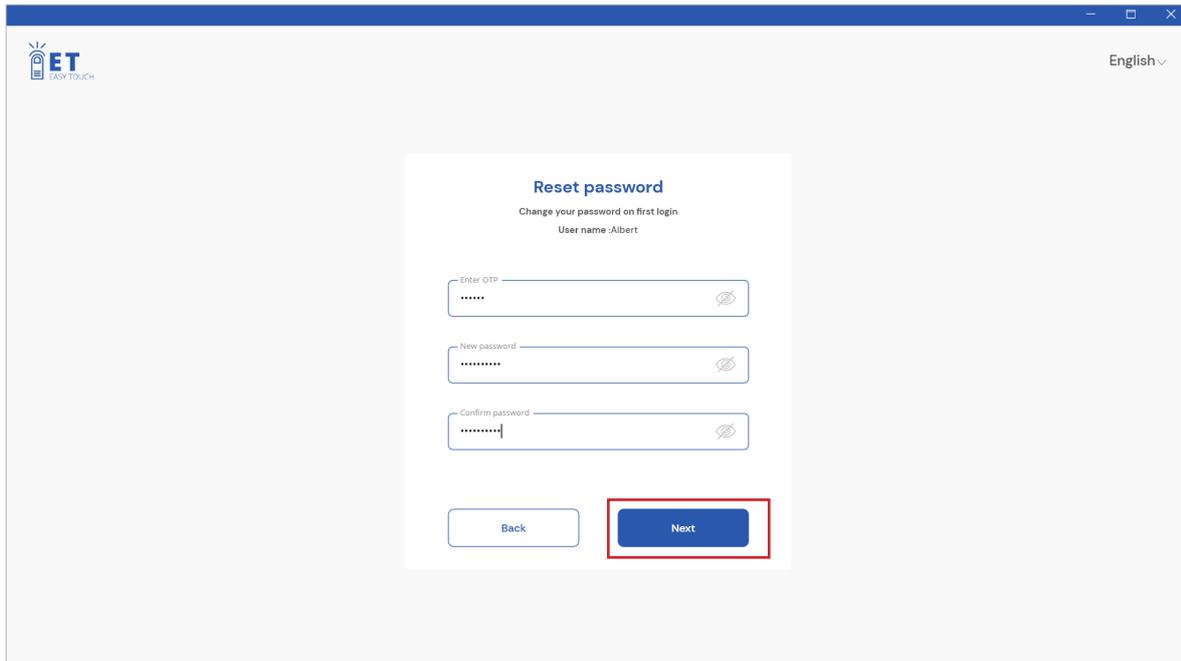


- The activated functions will available for usage.


2.4 Purchase licenses

Note:


- For the first time user, you may be able to see the screen to purchase or activate the licenses for the functions after logging in with the registered KERN account credentials.

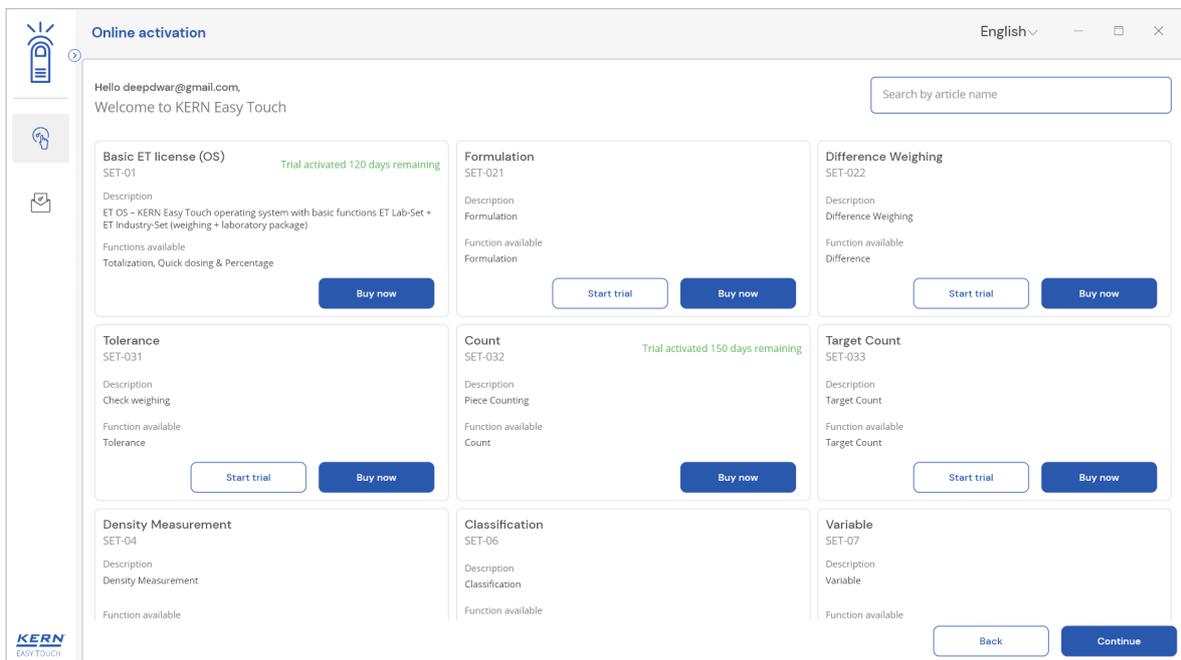

- For a non first time user (who have already activate the trial or purchased the licenses), the screen to purchase or activate the licenses for the functions would be available after logging in with admin account and navigating to settings> manage license> modify license

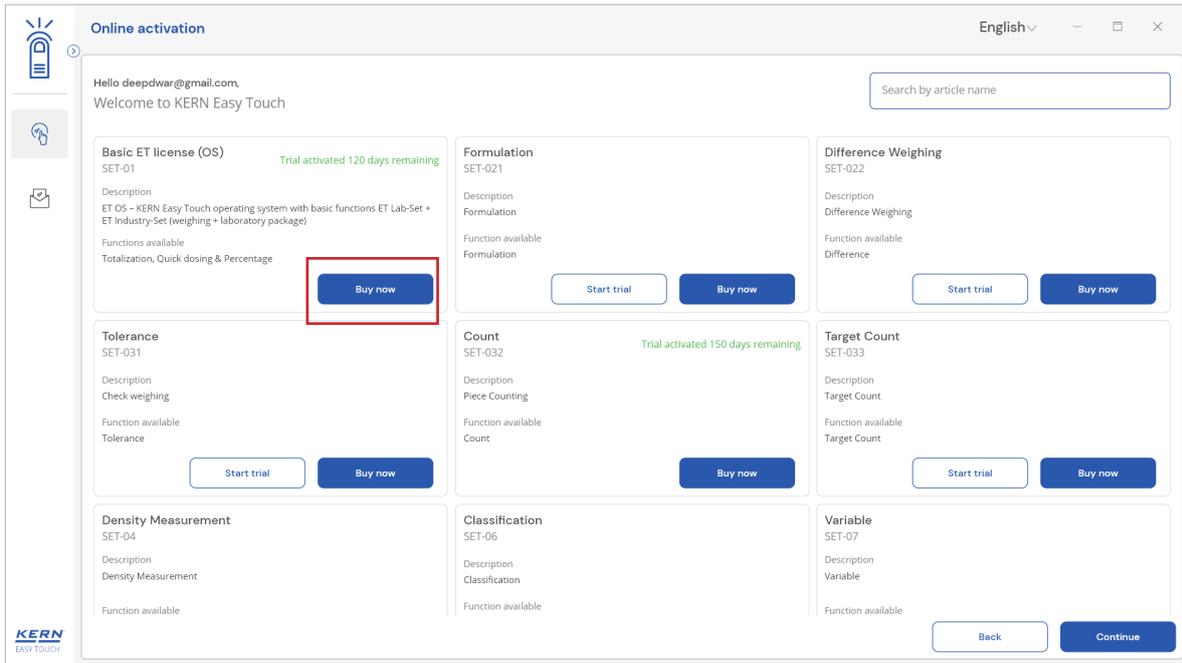


- Clicking on the “buy now” option will take you to the licensing website to purchase the required functions. (Please login with your KERN account credentials)

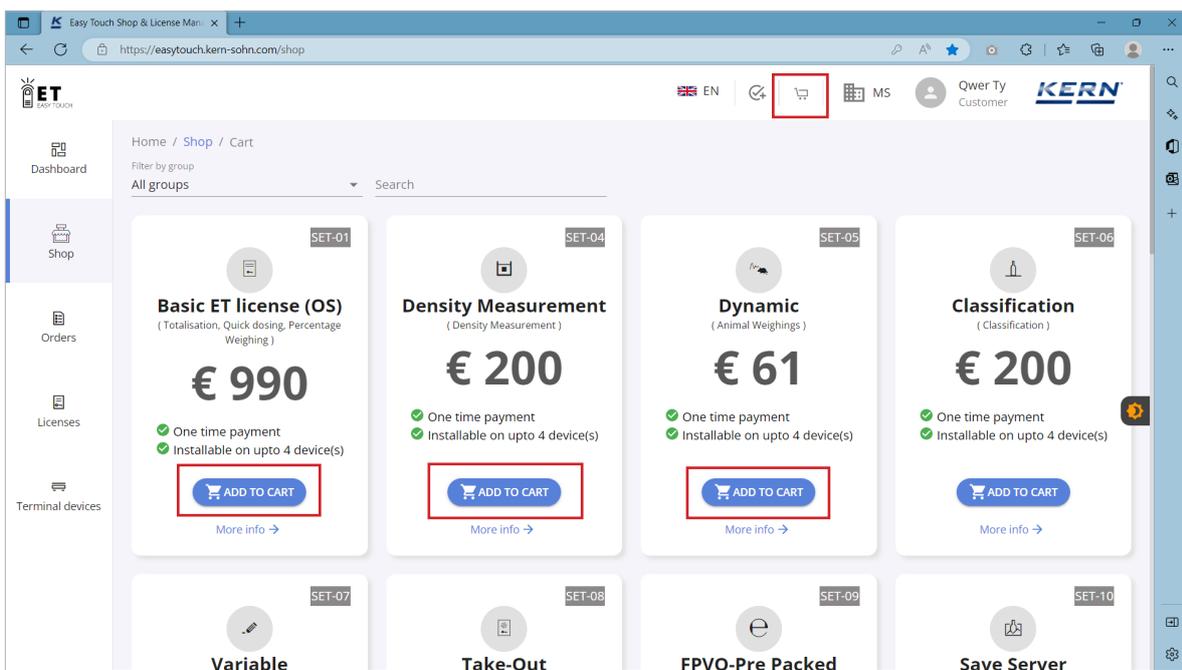


- Purchase your required functions by adding the licenses to cart and login to Easy Touch application to activate the required licenses.

2.5 License activation


- Once the license is purchased, you can now login the Easy Touch 2.0 application for activation.

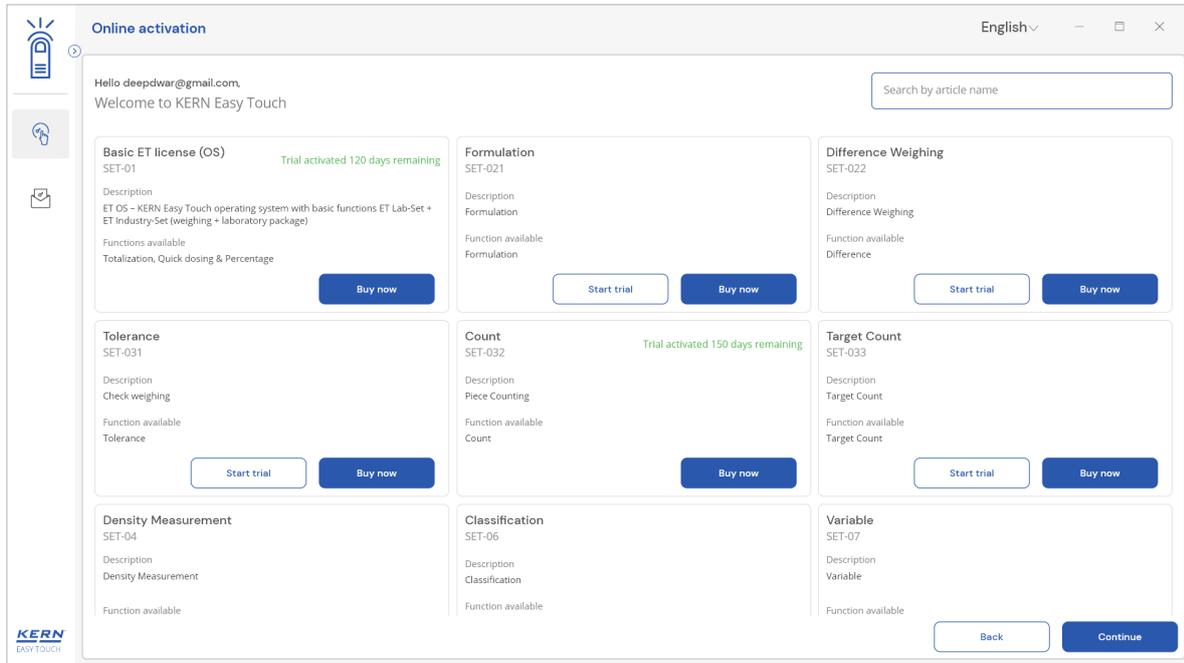


- The screen to purchase or activate the licenses for the functions would be available after logging in with admin account and navigating to settings> manage license> modify license. Enter the kern account credentials and activate the license.

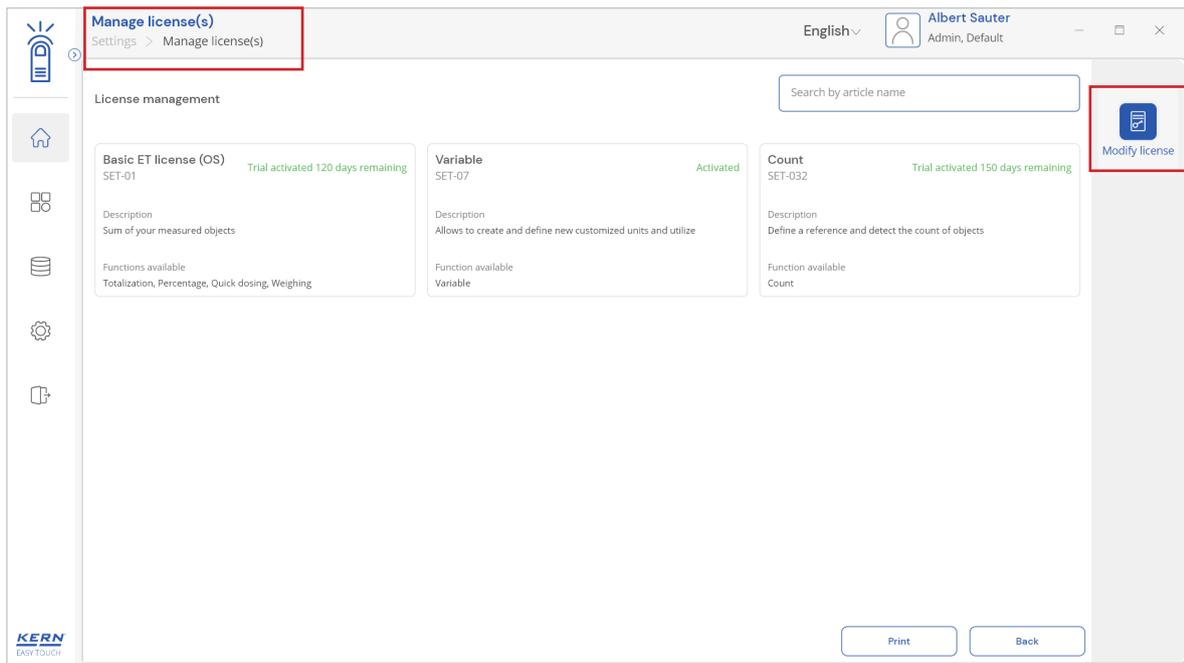

- Enter the created kern account username and password and click on login.

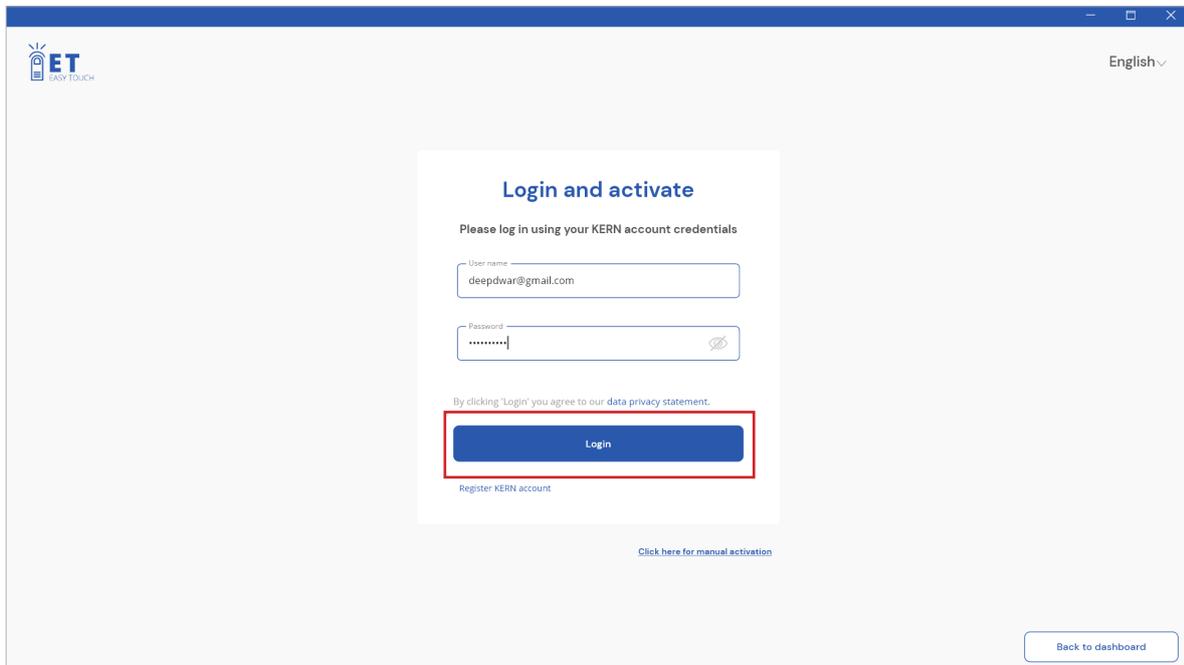


- When you click on the “login” button, you will be logged in to the easy touch application and redirected to the screen where you can activate your purchased license
- Click on “activate” for the license which you want to activate for the machine (SET - 01 is a mandatory function, to activate the other licenses, the basic os must be activated first.)

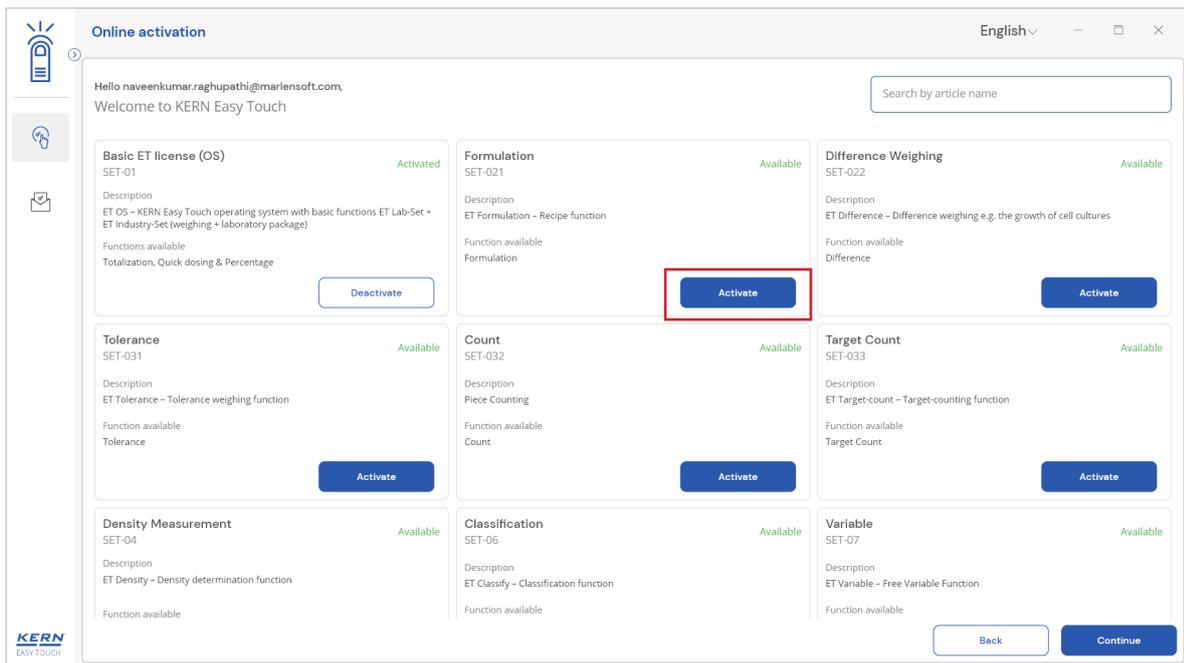


- Once the required licenses are activated in the machine click on “continue” to proceed.

English

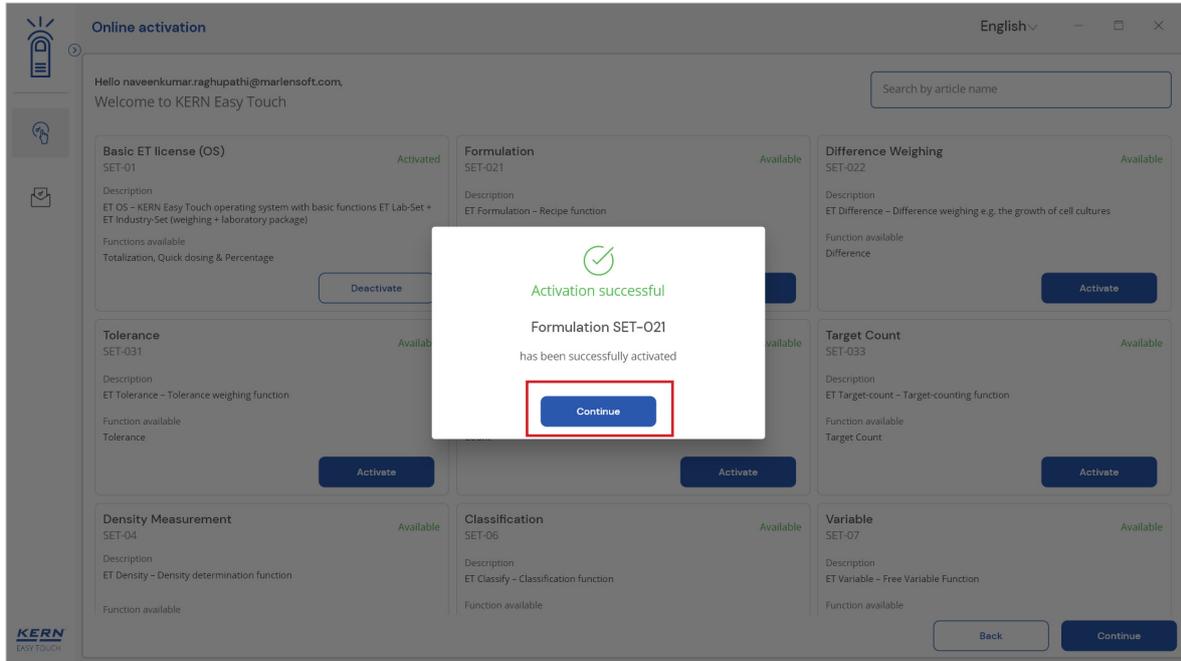

- If the license is activated then the status will be changed to “activated” and you will get the option to deactivate.

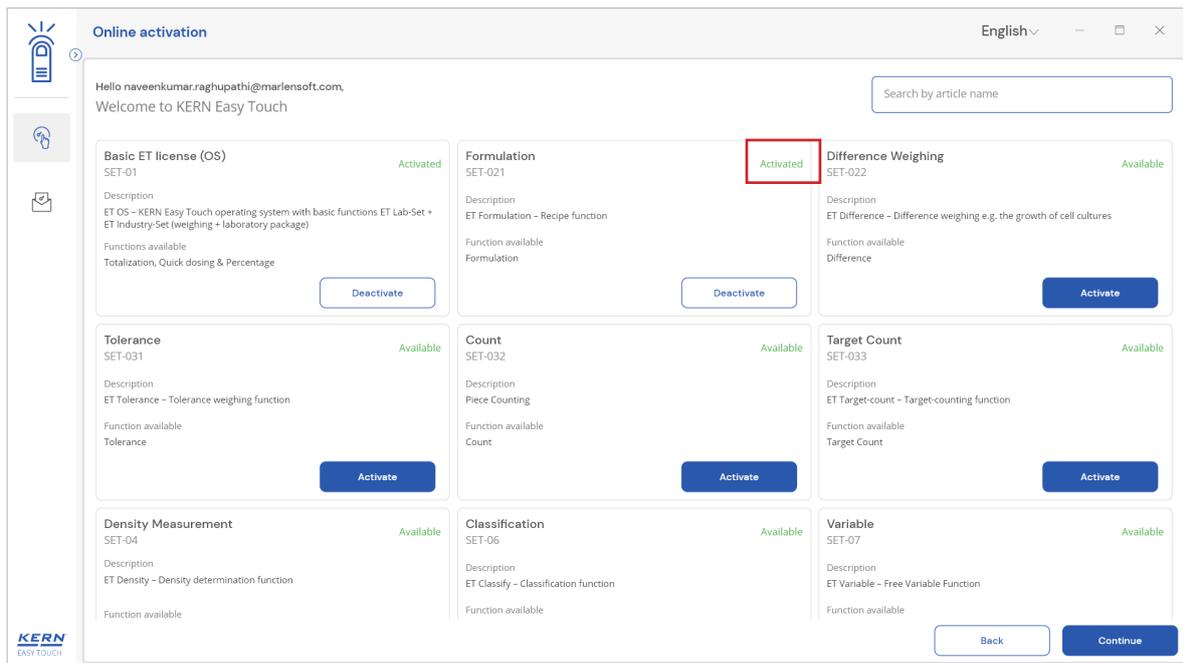

- Click on continue to utilise the functions

Online activation English

Hello naveenkumar.raghupathi@marlensoft.com
Welcome to KERN Easy Touch

Search by article name

Module Name	Article ID	Status	Action
Basic ET license (OS)	SET-01	Activated	Deactivate
Formulation	SET-021	Activated	Deactivate
Difference Weighing	SET-022	Available	Activate
Tolerance	SET-031	Available	Activate
Count	SET-032	Available	Activate
Target Count	SET-033	Available	Activate
Density Measurement	SET-04	Available	Activate
Classification	SET-06	Available	Activate
Variable	SET-07	Available	Activate

Back Continue

- Enter the admin credentials and login

Welcome to Easy Touch English


Enter your username and password to continue

User name: Albert

Password:


Login

English

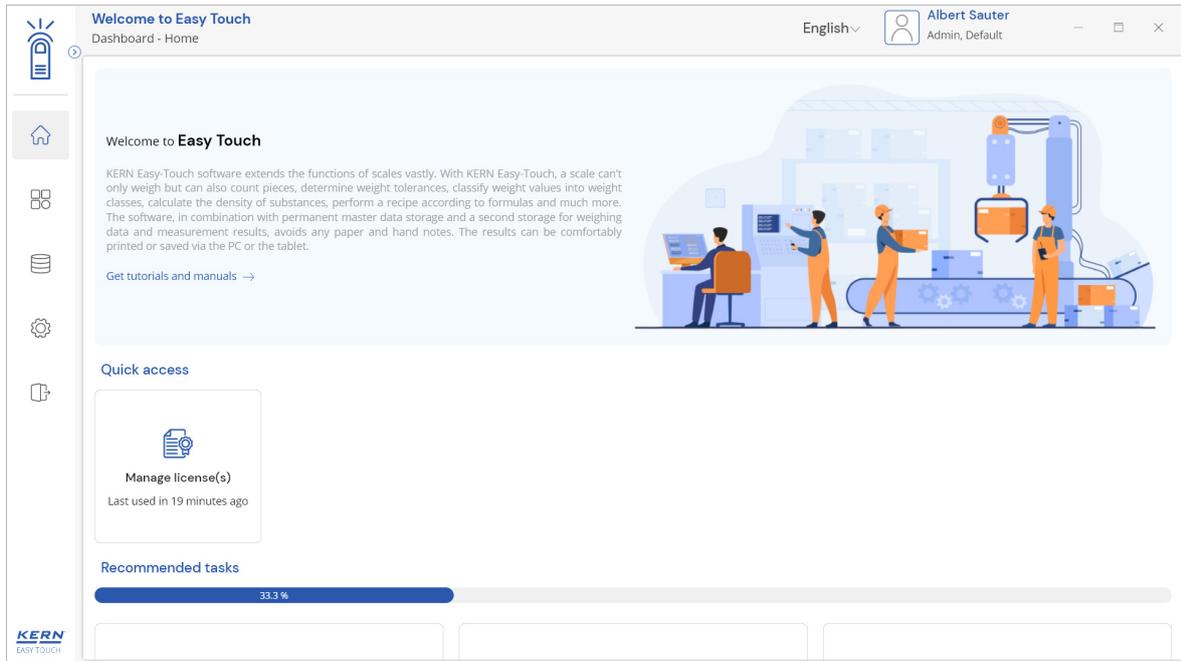

3.0 User with the registered Kern account

If the user has already a Kern account, then using the Kern account credentials the user can login directly once the application is installed and open

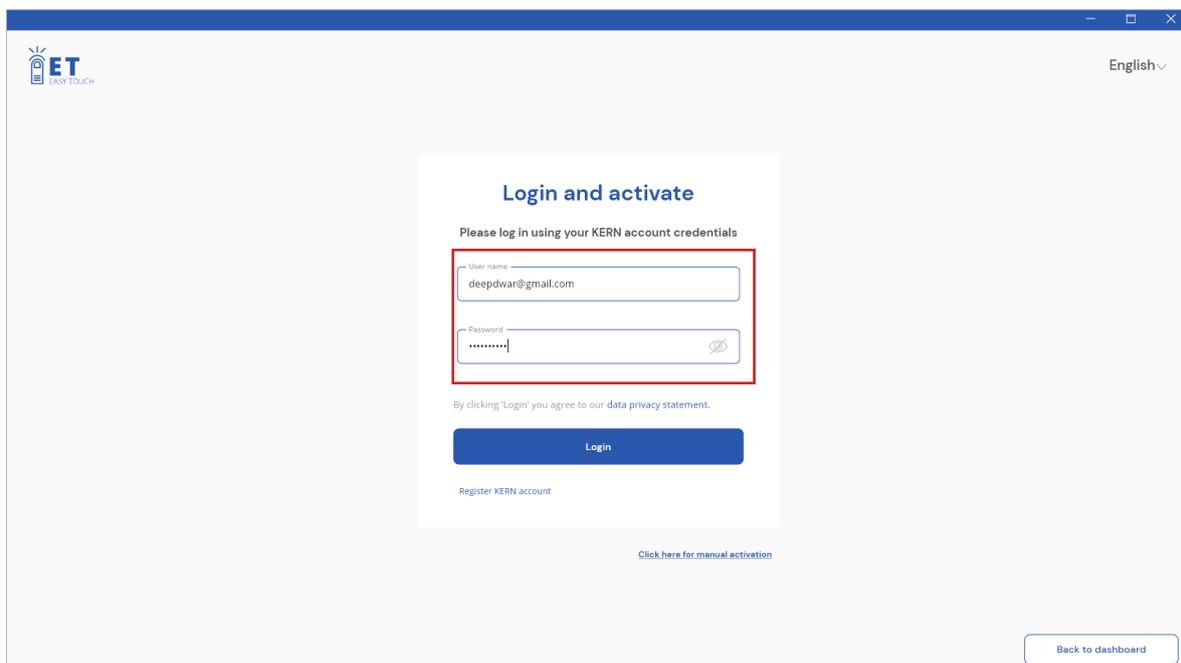

The user can activate the purchased licenses or purchase or activate the trial license or activate the purchased licenses by logging in with the KERN account credentials.

The following steps can be performed for using the Easy Touch functions according to the requirement,

- [Enable the trial license \(or purchase\)](#)
- [Admin creation for your system](#)
- [Purchase the functions](#)
- [Activation of purchased licenses](#)

Kindly refer the section 2.2, 2.3, 2.4 and 2.5 for the above steps.

The end