



**KERN & Sohn GmbH**

Ziegelei 1  
D-72336 Balingen  
E-mail: [info@kern-sohn.com](mailto:info@kern-sohn.com)

Tel.: +49-[0]7433-9933-0  
Fax: +49-[0]7433-9933-149  
Internet: [www.kern-sohn.com](http://www.kern-sohn.com)

# Installation manual

## Pedal

### KERN YKF-01

Version 1.2  
2018-08  
GB



YKF-01-IA-e-1812



# KERN YKF-01

Version 1.2 2018-08

## Installation manual Pedal

---

---



Only for KERN PCD and KERN PCB.

## 1 Installation

Attach the pedal to the device, as shown on the picture below.



## 2 Supplied elements

- Pedal
- Connector box
- 1 cable

## 3 Usage

The pedal is used to send weight values to a printer or a computer. Besides, it can be used to balance the scale.

### 3.1 Taring the scale

- ⇒ Turn the scale on using the **ON/OFF** button.
- ⇒ In weighing mode press and hold the **PRINT** button until the "**Unit**" menu is displayed.
- ⇒ Press the **MODE** button until the "**FOOt S**" parameter is displayed.
- ⇒ Confirm using the **SET** button, which will display the current setting.
- ⇒ Using the **Mode** button select the "**tAr**" setting and confirm it, using the **SET** button.
- ⇒ Place the scale container.
- ⇒ Press the pedal – the scale will display a zero result.

### 3.2 Printing or connecting to a computer



An appropriate cable compatible with the scale is required to transfer data. It should be connected between the connector box of the pedal and the printer or the computer.

- ⇒ Turn the scale on using the **ON/OFF** button.
- ⇒ In weighing mode press and hold the **PRINT** button until the "**Unit**" menu is displayed.
- ⇒ Press the **MODE** button until the "**FOOt S**" parameter is displayed.
- ⇒ Confirm using the **SET** button, which will display the current setting.
- ⇒ Using the **Mode** button select the "**Pr**" setting and confirm it, using the **SET** button.
- ⇒ Place the weighed material.
- ⇒ Press the pedal and the weighed value will be printed.