



KERN & Sohn GmbH

Ziegelei 1
D-72336 Balingen
E-Mail: info@kern-sohn.com

Phone: +49-[0]7433- 9933-0
Fax: +49-[0]7433-9933-149
Internet: www.kern-sohn.com

Installation instructions Battery holder

KERN PWS-A01

Version 1.0
2022-09
GB



TPWS-A01-A-IA-e-2210



KERN PWS-A01

Version 1.0 2022-09


Installation instructions Battery holder

1 Scope of delivery

- Battery holder
- Installation Instructions

2 Assembly

1. Loosen the two screws on the back of the balance and remove the cover

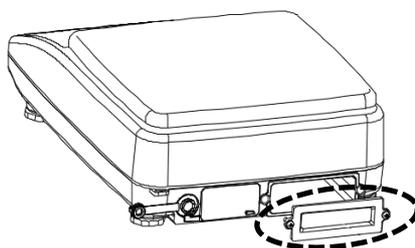


2. Insert the batteries / rechargeable batteries into the battery holder:
4 x 1.5 V AA

Pay attention to the polarity when inserting them



3. Slide the battery holder into the opening and tighten with the rotary knobs

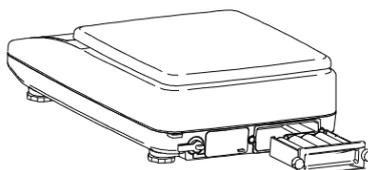

Please note that the rechargeable batteries cannot be charged using the balance.

3 Display of the charge level

Symbol	Description
	Charge level: normal
	Charge level: low
	Charge level: empty (flashing symbol)